LABCONCO CORPORATION

8811 Prospect Avenue
Kansas City, MO 64132
(800) 821-5525, (816) 333-8811
(816) 363-0130 fax
labconco@labconco.com

User's Manual

Protector[®] XL[™] Floor-Mounted (Walk-In) Laboratory Fume Hoods

Models

1124 Series, 4' Vertical
1125 Series, 5' Vertical
1126 Series, 6' Vertical
1128 Series, 8' Vertical & Horizontal
1120 Series, 10' Horizontal
1121 Series, 12' Horizontal
1122 Series, 16' Horizontal

To receive important product updates,
complete your product registration card
online at register.labconco.com

Please read the User's Manual before operating the equipment.

Copyright © 2021 Labconco Corporation. The information contained in this manual and the accompanying products are copyrighted and all rights reserved by Labconco Corporation. Labconco Corporation reserves the right to make periodic design changes without obligation to notify any person or entity of such change.

Warranty

Labconco Corporation provides a warranty to the original buyer for the repair or replacement of parts and reasonable labor as a result of normal and proper use of the equipment with compatible chemicals. Broken glassware and maintenance items, such as filters, gaskets, light bulbs, finishes and lubrication are not warranted. Excluded from warranty are products with improper installation, erratic electrical or utility supply, unauthorized repair and products used with incompatible chemicals.

The warranty for Protector® XL™ Floor-Mounted (Walk-In) Laboratory Fume Hoods will expire one year from date of installation or two years from date of shipment from Labconco, whichever is sooner. Warranty is non-transferable and only applies to the owner (organization) of record.

Buyer is exclusively responsible for the set-up, installation, verification, decontamination or calibration of equipment. This limited warranty covers parts and labor, but not transportation and insurance charges. If the failure is determined to be covered under this warranty, the dealer or Labconco Corporation will authorize repair or replacement of all defective parts to restore the unit to operation. Repairs may be completed by 3rd party service agents approved by Labconco Corporation. Labconco Corporation reserves the rights to limit this warranty based on a service agent's travel, working hours, the site's entry restrictions and unobstructed access to serviceable components of the product.

Under no circumstances shall Labconco Corporation be liable for indirect, consequential, or special damages of any kind. This warranty is exclusive and in lieu of all other warranties whether oral, or implied.

Returned or Damaged Goods

Do not return goods without the prior authorization from Labconco. Unauthorized returns will not be accepted. If your shipment was damaged in transit, you must file a claim directly with the freight carrier. Labconco Corporation and its dealers are not responsible for shipping damages.

The United States Interstate Commerce Commission rules require that claims be filed with the delivery carrier within fifteen (15) days of delivery.

Limitation of Liability

The disposal and/or emission of substances used in connection with this equipment may be governed by various federal, state, or local regulations. All users of this equipment are required to become familiar with any regulations that apply in the user's area concerning the dumping of waste materials in or upon water, land, or air and to comply with such regulations. Labconco Corporation is held harmless with respect to user's compliance with such regulations.

Contacting Labconco Corporation

If you have questions that are not addressed in this manual, or if you need technical assistance, contact Labconco's Customer Service Department or Labconco's Product Service Department at 1-800-821-5525 or 1-816-333-8811, between the hours of 7:30 a.m. and 5:30 p.m., Central Standard Time.

TABLE OF CONTENTS

CHAPTER 1: INTRODUCTION	1
About This Manual	3
Typographical Conventions (Conventions Typographiques)	4
Your Next Step	5
CHAPTER 2: PREREQUISITES	7
Location Requirements	8
Support Requirements	8
Exhaust Requirements (Exigences électriques)	8
Electrical Requirements	10
Service Line Requirements	10
Space Requirements	10
CHAPTER 3: GETTING STARTED	11
Unpacking Your Floor-Mounted Fume Hood	12
Removing the Shipping Skid	12
Sash Weight Release	13
Disassembly & Reassembly of the XL Floor-Mounted Hood	13
Lower Base Installation of the XL Floor-Mounted Hood	13
Upper Cabinet Installation of the XL Floor-Mounted Hood	14
Vertical-Rising Sash Installation	15
Lower Sash Track & Threshold Installation	19
Horizontal-Sliding Sash Installation	20
Corner Post Installation for the Floor-Mounted Hoods	21
Baffle Installation for the Floor-Mounted Hoods	21
Header Installation for the Floor-Mounted Hoods	23
Front Panel Installation for the Floor-Mounted Hoods	23
Connecting to the Hood Exhaust System	24
Connecting the Electrical Supply Source	24
Connecting the Service Lines to the Protector Hood	27
Sealing the Protector Hood to the Floor	28
Certifying the Protector Fume Hood	28


CHAPTER 4: PERFORMANCE FEATURES AND SAFETY PRECAUTIONS	29
Performance Features	29
Safety Precautions (Consignes de sécurité)	32
CHAPTER 5: USING YOUR PROTECTOR XL	36
Operating the Vertical-Rising Sash	36
Operating the Horizontal-Sliding Sashes	36
Operating the A-Style Combination Sash	36
Operating the Blower	37
Operating the Lights	37
Working in your Protector Fume Hood	37
CHAPTER 6: MAINTAINING YOUR PROTECTOR XL	39
Service Safety Precautions (Mesures de sécurité des services)	39
Routine Maintenance Schedule	40
Routine Service Operations	41
CHAPTER 7: MODIFYING YOUR PROTECTOR XL	42
Installing Work Surfaces	42
Installing Ceiling Enclosures Above the Fume Hood	42
Installing Additional Service Fixtures	45
Installing Guardian Airflow Monitors	46
Floor Mount Distillation Grids – Field Installation	46
Sash Stop Kit – Field Installation	47
Installing an Electrical Duplex Outlet (Installation d'une sortie Duplex électrique)	47
8' Dual Exhaust Adapter	47
CHAPTER 8: TROUBLESHOOTING	48
APPENDIX A: PROTECTOR XL COMPONENTS	51
APPENDIX B: PROTECTOR XL DIMENSIONS	55
APPENDIX C: PROTECTOR XL SPECIFICATIONS (ANNEXE C: SPÉCIFICATIONS PROTECTEUR DU PREMIER MINISTRE)	57
APPENDIX D: SERIAL NUMBER TAG DESCRIPTION (ANNEXE D: DESCRIPTION DE SÉRIE TAG NUMBER)	60
APPENDIX E: REFERENCES	66
APPENDIX F: DISASSEMBLY AND REASSEMBLY INSTRUCTIONS	69
APPENDIX G: VAV FLOOR-MOUNTED OPEN AREA GUIDELINES	72

CHAPTER 1


INTRODUCTION

Congratulations on your purchase of a Labconco Protector® XL™ Floor-Mounted (Walk-In) Laboratory Fume Hood. Your Protector Laboratory Floor-Mount Fume Hood is designed to protect the laboratory scientist who must transport equipment and products for experimentation and testing under a laboratory hood. It is the result of Labconco's more than 60 years of experience in manufacturing fume hoods, and users like you suggested many of its features to us. The Vertical Floor-Mounted Hood is rated high performance low flow at 60 fpm and the Horizontal Floor-Mounted Hood is listed at 80 fpm.

The Labconco Protector XL Floor-Mounted Fume Hood has been engineered to provide maximum visibility in a laboratory, and effectively contain toxic, noxious, or other harmful materials when properly installed. The Protector XL offers many unique features to enhance safety, performance, and visibility. To take full advantage of them, please acquaint yourself with this manual and keep it handy for future reference. If you are unfamiliar with how floor-mount fume hoods operate, please review *Chapter 4: Performance Features and Safety Precautions* before you begin working in the fume hood. Even if you are an experienced fume hood user, please review *Chapter 5: Using Your Protector XL*, which describes your Protector Floor-Mounted Hood's features so that you can use the hood efficiently.

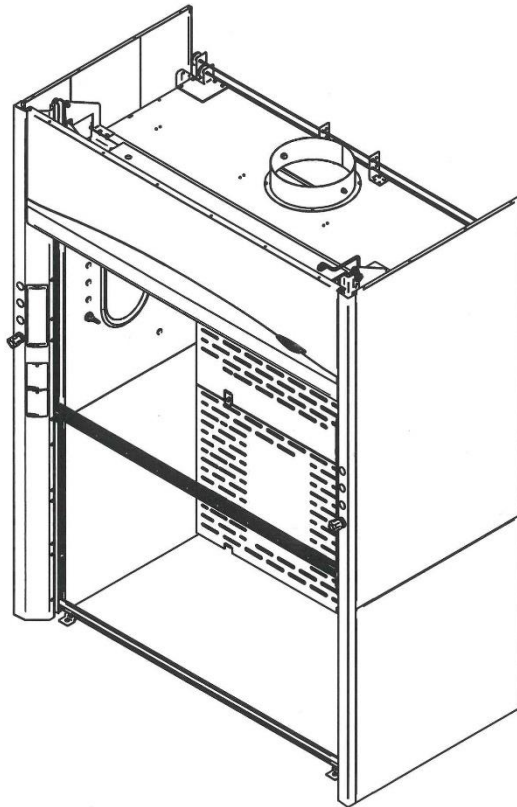**Figure 1-1
Vertical Floor-Mount**

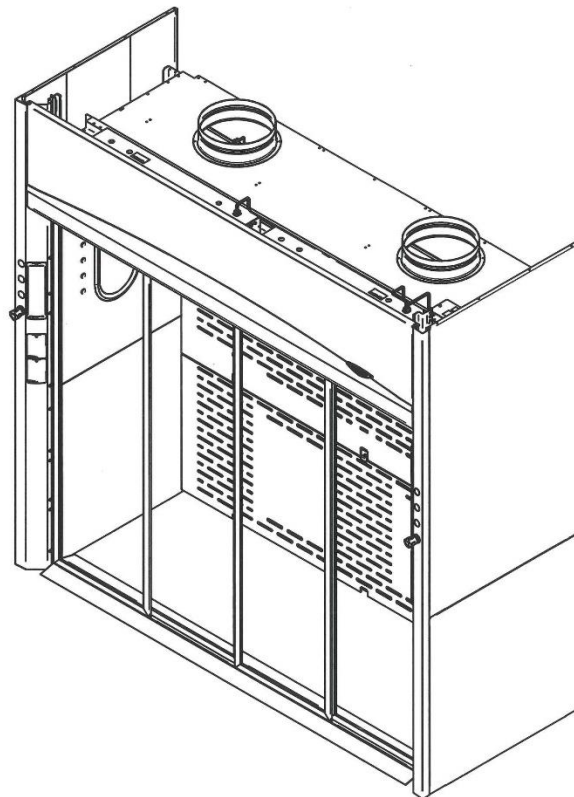**Figure 1-2
Horizontal Floor-Mount**

About This Manual

This manual is designed to help you learn how to install, use, and maintain your laboratory fume hood. Instructions for installing optional equipment on your hood are also included.

Chapter 1: Introduction provides a brief overview of the floor-mounted laboratory fume hood, explains the organization of the manual, and defines the typographical conventions used in the manual.

Chapter 2: Prerequisites explains what you need to do to prepare your site before you install your floor-mounted laboratory fume hood. Electrical and service requirements are discussed.

Chapter 3: Getting Started contains the information you need to properly unpack, inspect, install, and certify your floor-mounted laboratory fume hood.

Chapter 4: Performance Features and Safety Precautions explains how the Protector operates and the appropriate precautions you should take when using the floor-mounted fume hood.

Chapter 5: Using Your Protector XL Floor-Mounted Hood discusses the basic operation of your fume hood. Information on how to prepare, use and shut down your Protector Hood are included.

Chapter 6: Maintaining Your Protector XL Floor-Mounted Hood explains how to perform routine maintenance on your fume hood.

Chapter 7: Modifying Your Protector XL Floor-Mounted Hood explains how to modify the fume hood or add accessories.

Chapter 8: Troubleshooting contains a table of problems you may encounter while using your laboratory fume hood including the probable causes of the problems and suggested corrective actions.

Appendix A: Protector XL Floor-Mounted Hood Components contains labeled diagrams of all of the components of the fume hoods.

Appendix B: Protector XL Floor-Mounted Hood Dimensions contains comprehensive diagrams showing all of the dimensions for the laboratory fume hoods.

Appendix C: Protector XL Floor-Mounted Hood Specifications contains the electrical requirements for laboratory fume hood. Wiring diagrams are also included.

Appendix D: Serial Number Tag Description provides current rating code used on serial number tag.

Chapter 1: Introduction

Appendix E: References lists the various resources available that deal with laboratory fume hoods.

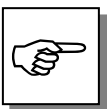
Appendix F: Disassembly and Reassembly Instructions lists the required tools as well as step by step instructions for disassembling and reassembling the Protector XL Hood.

Appendix G: VAV Floor-Mounted Open Area Guidelines to properly assess fume hood effective area whenever variable air volume (VAV) mechanical systems are installed.

Typographical Conventions

Recognizing the following typographical conventions will help you understand and use this manual:

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- 

 - A number icon precedes information that is specific to a particular model of laboratory fume hood. The 4' icon indicates the text is specific to the 4-foot wide model. The 5' icon indicates the text is specific to the 5-foot model, etc.
- 

- 
 - The S icon indicates the text is specific to the standard model.
- 
 - The A icon indicates the text is specific to the A-Style Combination Sash Model.
- 
 - CAUTION – See Manual. When this symbol is on a fume hood it indicates a caution that is detailed in this manual.
 - PRUDENCE – Consulter le Manuel. Quand ce symbole est sur une hotte aspirante, il indique une prudence qui est détaillée dans ce manuel.
- 
 - CAUTION – Hot Surface.
 - AVERTIR – Surface Chaude
- 
 - CAUTION – See Manual. This symbol on the fume hood indicates the possibility of a pinch hazard.
 - PRUDENCE – Consulter le Manuel. Ce symbole sur la hotte indique la possibilité d'un risque de pincement.

Your Next Step

If your Fume Hood needs to be installed, proceed to *Chapter 2: Prerequisites* to ensure your installation site meets all of the requirements. Then, go to *Chapter 3: Getting Started* for instructions on how to install your laboratory fume hood and make all of the necessary connections.

If you would like to review how laboratory fume hoods operate, go to *Chapter 4: Performance Features and Safety Precautions*.

For information on the operational characteristics of your laboratory fume hood, go to *Chapter 5: Using Your Protector XL Floor-Mounted Hood*.

If your laboratory fume hood is installed and you need to perform routine maintenance on the cabinet, proceed to *Chapter 6: Maintaining Your Protector XL Floor-Mounted Hood*.

Chapter 1: Introduction

For information on making modifications to the configuration of your fume hood, go to *Chapter 7: Modifying Your Protector XL Floor-Mounted Hood*.

Refer to *Chapter 8: Troubleshooting* if you are experiencing problems with your fume hood.

CHAPTER 2

PREREQUISITES

Before you install your floor-mount laboratory fume hood, you need to prepare your site for installation. Carefully examine the location and floor where you intend to install your floor-mounted hood. You must be certain that the area is level and of solid construction. In addition, a dedicated source of electrical power must be located near the installation site.

Carefully read this chapter to learn the requirements for your installation site:

- The location requirements.
- The support requirements.
- The exhaust requirements.
- The electrical power requirements.
- The service line requirements.
- The space requirements.

Refer to *Appendix B: Protector XL Floor-Mounted Hood Dimensions* for complete fume hood dimensions.

Refer to *Appendix C: Protector XL Floor-Mounted Hood Specifications* for complete laboratory fume hood electrical and environmental conditions, specifications and requirements.

Location Requirements

The floor-mounted fume hood should be located away from traffic patterns, doors, windows, fans, ventilation registers, and any other air-handling device that could disrupt its airflow patterns. All windows in the room should be closed.

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Support Requirements

DO NOT install the fume hood on a cart, dolly, or mobile bench. **ALL** Protector Hood installations must be permanent and stationary. The Protector XL Floor-Mounted Hood is usually mounted directly on room floor surfaces, or an optional chemical resistant work surface.

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Exhaust Requirements

The exhaust duct connection has been designed for 12" nominal duct (12.75" OD) to allow for minimum static pressure loss while operating at 60 to 100 fpm face velocities. The 12" diameter exhaust duct also allows for proper transport velocities away from the hood in the 1000 fpm to 2500 fpm range. The proper exhaust volume and static pressure loss are listed next for each hood model:

Projected Airflows and Static Pressure

Hood Type	Face	Velocity (fpm)	Airflow Volumetric Rate (CFM) @ Static Pressure (inches of water)			
	Top Sash Full open at 64"	Top Sash open 62.5% at 54"	4' Hood	5' Hood	6' Hood	8' Hood
Vertical						
Floor						
Mount						
Walk-In	100	160	725, 0.27"	955, 0.34"	1180, 0.46"	1640, 0.31"
XL	80	128	580, 0.17"	765, 0.22"	945, 0.29"	1310, 0.20"
W/Bypass	62.5	100	450, 0.11"	595, 0.13"	735, 0.18"	1025, 0.12"
Block	60	96	435, 0.10"	575, 0.12"	710, 0.17"	985, 0.11"
	50	80	365, 0.07"	480, 0.09"	590, 0.11"	820, 0.08"
	40	64	290, 0.05"	380, 0.06"	470, 0.08"	655, 0.05"
	38	60	270, 0.04"	360, 0.05"	440, 0.07"	615, 0.04"

Note: Bottom Sash is closed and only opened for loading.

Projected Airflows and Static Pressure

Hood Type	Face	Velocity (fpm)	Doors Open	2 Doors	1 Door	1 Door	1 Door
			Airflow Volumetric Rate (CFM) @ Static Pressure (inches of water)				
Horizontal	Horizontal			8' Hood	10' Hood	12' Hood	16' Hood
Floor	Sash at						
Mount	Full Open						
Walk-In							
XL	100		2150, 0.53"	1450, 0.22"	1765, 0.30"	2400, 0.18"	
W/Bypass	80		1720, 0.34"	1160, 0.14"	1410, 0.19"	1920, 0.12"	
Block							
			2 Doors Open	1 Door Open	1 Door Open	1 Door Open	

Proper blower selection for vertical and horizontal floor-mounted hoods can be determined from these exhaust requirements and the total system static pressure loss. Contact Labconco Customer Service for assistance in sizing a blower system. Refer to *Appendix G: VAV Floor-Mounted Open Area Guidelines* to properly assess the fume hood area opening, allowing the variable air volume (VAV) mechanical system to be programmed at various sash positions.

Electrical Requirements

The Protector XL Floor-Mounted Hood models feature internal wiring for the LED light assembly and light switch. All internal wiring is terminated at the single point wiring junction box for hook-up by a qualified electrician. The blower switch, and light switch wires are also terminated at the single point wiring junction box for hook-up by a qualified electrician. Refer to *Chapter 3: Getting Started* and *Appendix C: Protector XL Specifications* for the wiring diagram for proper electrical installation.

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Service Line Requirements

All service lines to the laboratory fume hood should be ¼ inch outside diameter, copper (brass for natural gas), and equipped with an easily accessible shut-off valve, should disconnection be required. Recommended operating pressure is 40 PSI, with a maximum allowable pressure of 200 PSI. Consider a pressure regulator to reduce line pressure to 40 PSI. Please check with local codes for other requirements.

Space Requirements

The dimensions for the different models are shown in *Appendix B: Protector XL Floor-Mounted Hood Dimensions*.

CHAPTER 3

GETTING STARTED

Now that the site for your floor-mounted fume hood is properly prepared, you are ready to unpack, inspect, install, and certify your unit. Read this chapter to learn how to:

- Unpack and move your floor-mounted hood.
- Set up the floor-mounted fume hood on the floor.
- Connect to an exhaust system.
- Connect the electrical supply source.
- Connect the service lines.
- Sealing the Protector Hood to the floor.
- Arrange certification of your Protector Hood.

Depending upon which model you are installing, you may need common plumbing and electrical installation tools in addition to 5/16", 3/8", 7/16", and 1/2" wrenches, ratchets, sockets, a nut driver set, a flat-blade screwdriver, a Phillips screwdriver, and a carpenter level to complete the instructions in the chapter.

The Protector XL Hood models weigh between 700 to 3500 lbs. (318-1593 kg). The multiple shipping pallets allow for lifting with a mechanical lift truck or floor jack. If you must lift the fume hood manually, follow safe-lifting guidelines.

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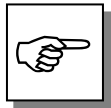
Unpacking Your Floor-Mounted Fume Hood


The United States Interstate Commerce Commission rules require that claims be filed with the delivery carrier within fifteen (15) days of delivery.

Your Protector XL Floor-Mounted (Walk-In) Hood has been shipped to you as ten main component assemblies located on three shipping skids. The ten main component assemblies consist of the lower base, upper cabinet, sashes, lower sash track, corner posts, service fixtures, electrical connections, baffles, header, and front panel.

Carefully remove the shrink-wrap or carton on your fume hood and inspect it for damage that may have occurred in transit. If your unit is damaged, notify the delivery carrier immediately and retain the entire shipment intact for inspection by the carrier.

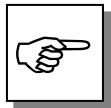DO NOT RETURN GOODS WITHOUT THE PRIOR AUTHORIZATION OF LABCONCO. UNAUTHORIZED RETURNS WILL NOT BE ACCEPTED.

IF YOUR HOOD WAS DAMAGED IN TRANSIT, YOU MUST FILE A CLAIM DIRECTLY WITH THE FREIGHT CARRIER. LABCONCO CORPORATION AND ITS DEALERS ARE NOT RESPONSIBLE FOR SHIPPING DAMAGES.

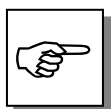
Do not discard the shipping skid or packing material for your floor-mounted fume hood until you have checked all of the components and installed and tested the unit. **Many components of the floor-mounted hood are shipped loose such as baffles and corner posts and do not discard. See Figure 1-1, Figure 1-2, Figure 3-1, Figure 3-7, Figure B-1 and Figure B-2 that shows proper XL Baffle installation.** Do not remove the fume hood from its shipping skids until it is ready to be placed into its final location. Move the unit by placing a flat, low dolly under the shipping skid, or by using a floor jack.

Do not move the hood by tilting it onto a hand truck.

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Removing the Shipping Skid


LEAVE THE FUME HOOD ATTACHED TO ITS SHIPPING SKID UNTIL IT IS AS CLOSE TO ITS FINAL LOCATION AS POSSIBLE. MOVE THE HOOD BY USING A SUITABLE FLOOR JACK, OR BY PLACING A FURNITURE DOLLY UNDERNEATH THE SKID. DO NOT MOVE THE HOOD BY TILTING IT ONTO A HAND TRUCK.

After you verify the fume hood components, move your hood to the location where you want to install it. Then, follow the steps listed next to remove the separate shipping skids from the lower base and upper cabinet.

1. Remove the side panels by unscrewing the Phillips screws.
2. Find the hardware (bolts, washers, nuts) that attach the fume hood to the skid and remove the hardware. Some hardware is on the sides and some is on the back.

Sash Weight Release

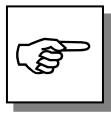On Floor-Mounted Hood models with vertical-rising sashes, the sash weights have been secured to the shipping skid. Remove the weights from the skid and attach them to the respective sash cables using the hooks provided.

NOTE: THE SASH WEIGHT ITSELF WAS INDIVIDUALLY MATCHED FOR THIS SPECIFIC HOOD AND SHOULD NOT BE EXCHANGED ON ANY OTHER UNIT.

Disassembly and Reassembly of the Protector XL Floor-Mounted Hood

There may be some disassembly and reassembly work, due to the large physical size of your laboratory hood and the ability to maneuver it into your laboratory. This is specific to each customer.

Lower Base Installation of the Protector XL Floor-Mounted Hood

V

The lower base features a panelized liner surrounded by an epoxy-coated steel and galvanized steel framework. Both the left and right side panels should be removed to allow for the correct placement of the lower cabinet assembly. Temporarily remove both vertical-rising sashes shipped with the lower base (on vertical-rising sash models only). Position the lower base on a level floor surface. Place the vertical-rising sashes back in place (on vertical-rising sash models only).

NOTE: If an optional floor surface is purchased, it should be positioned on the existing floor prior to placement of the lower cabinet installation. See Figure 3-1.

Upper Cabinet Installation of the Protector XL Floor-Mounted Hood

Remove the upper cabinet assembly side panels and front panels prior to cabinet placement. Then place the upper cabinet assembly on top of the lower base assembly, being careful to clear the lower base assembly during placement. To prevent spillage from seeping between the two liner assemblies, run a bead of white RTV sealant between the sections once they have been properly aligned. Mount the lower base to the upper cabinet using the 1/4-20 hex head screws, lockwashers, and nuts included in your hood package. See Figure 3-2.

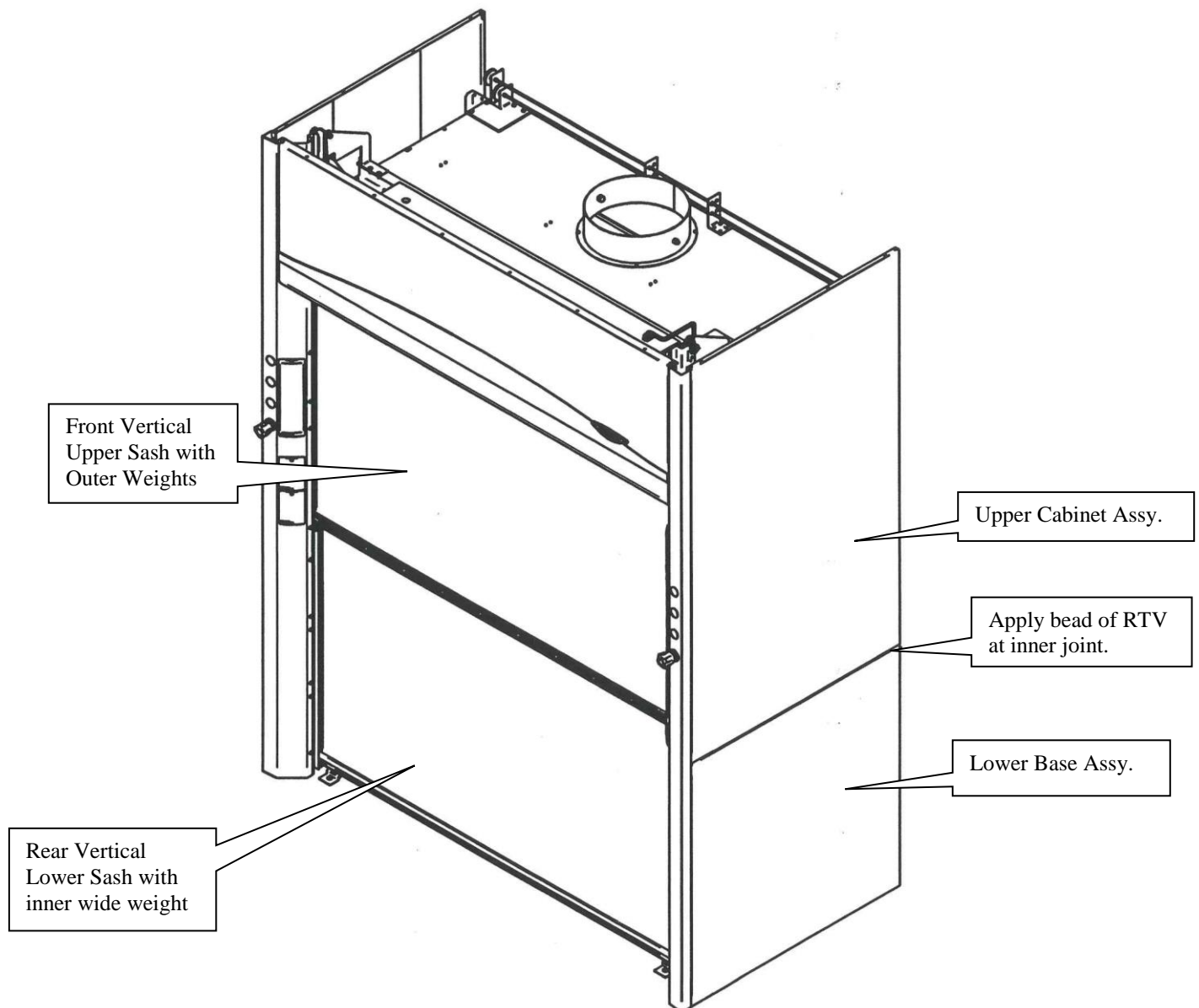
Figure 3-1

Vertical-Rising Sash Installation for the 4', 5', 6', & 8' Floor-Mounted Hoods

V

The Vertical-Rising Sash Floor-Mounted Hood has two vertical-rising sashes. The rear vertical-rising sash moves from the floor to full open and picks up the front vertical-rising sash, which travels from the midpoint to full open. The vertical rising sashes are shipped captured between the hood side frame assemblies of the lower base assembly. Both upper and lower sash tracks are attached to the side frames. The hood should be at a pre-assembled state with the top section bolted to the base, and the sashes still captured between the two side frames. Bring the sides back and adjust the sash tracks left and right to assure that each sash will work properly and tighten all hardware.

Now that the sashes are in place, the counterbalance weights must be installed. The large single sheet metal sash weight is installed in the middle and attached to the lower rear sash. The two individual sash weights counterbalance the front vertical-rising sash; these weights have rollers and ride in the rear sash tracks that straddle the middle sheet metal sash weight. Install the rear weight tracks for the individual sash weights with the #10-24 screws and #10-24 KEPS nuts supplied. See Figures 3-2, 3-3, and 3-4. It is vitally important that the front upper and lower sash tracks are fully aligned and the strap plates are used to secure this alignment. See Figure 3-3 and Figure 3-4. Once the sashes are aligned and operating freely, secure the lower base to the floor with sealant or fasteners. (Note: Hardware not supplied due to installation variables.) Securing the lower base will ensure the sashes will work freely.

As a last step, install the four rubber bumpers with #6 screws supplied. One set of rubber bumpers prevents the front upper sash from traveling too low. The second set of rubber bumpers prevent the rear lower sash from touching the floor and provides a good sweep of clean air across the floor. See Figure 3-3.

Chapter 3: Getting Started

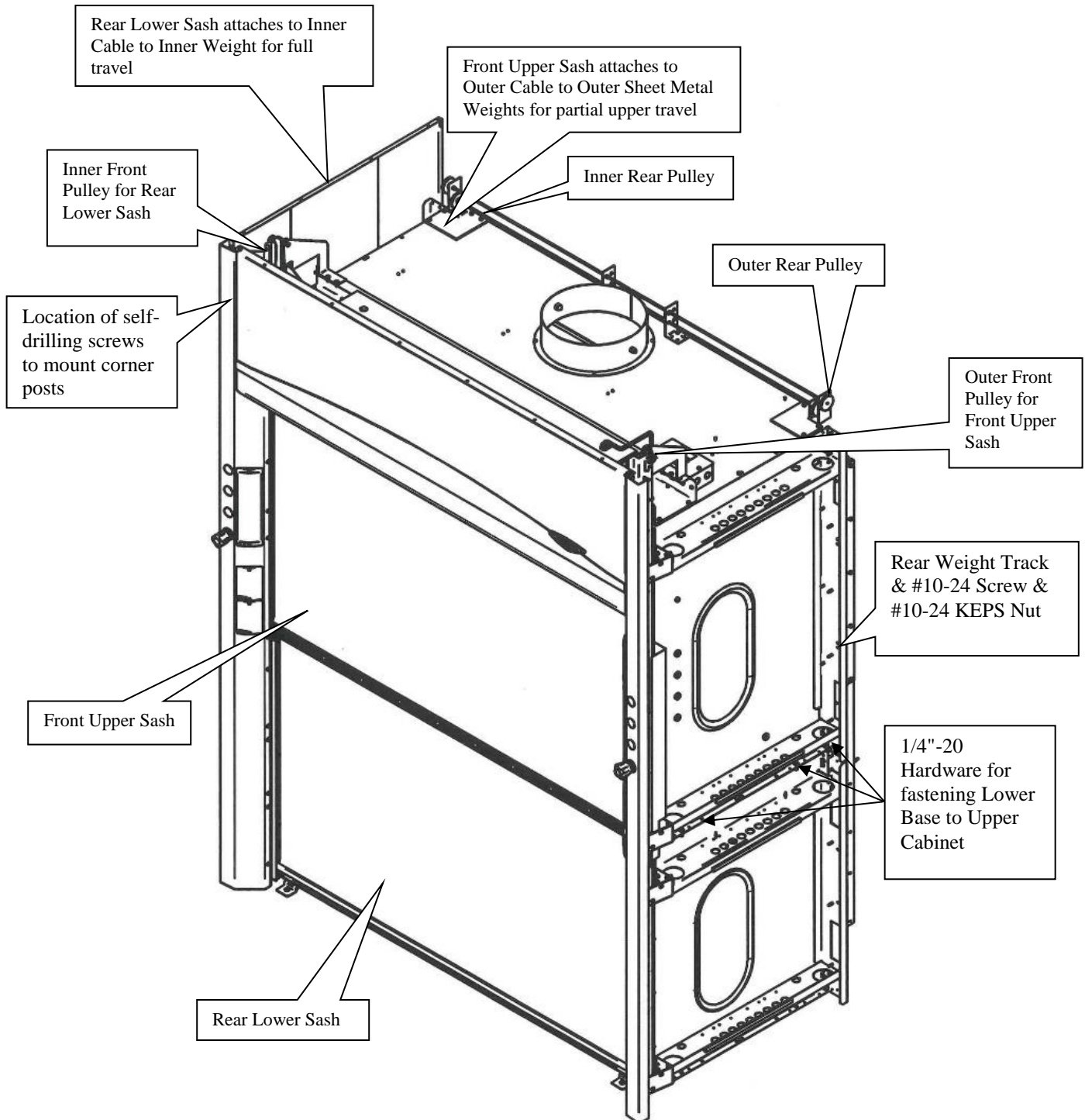
Figure 3-2


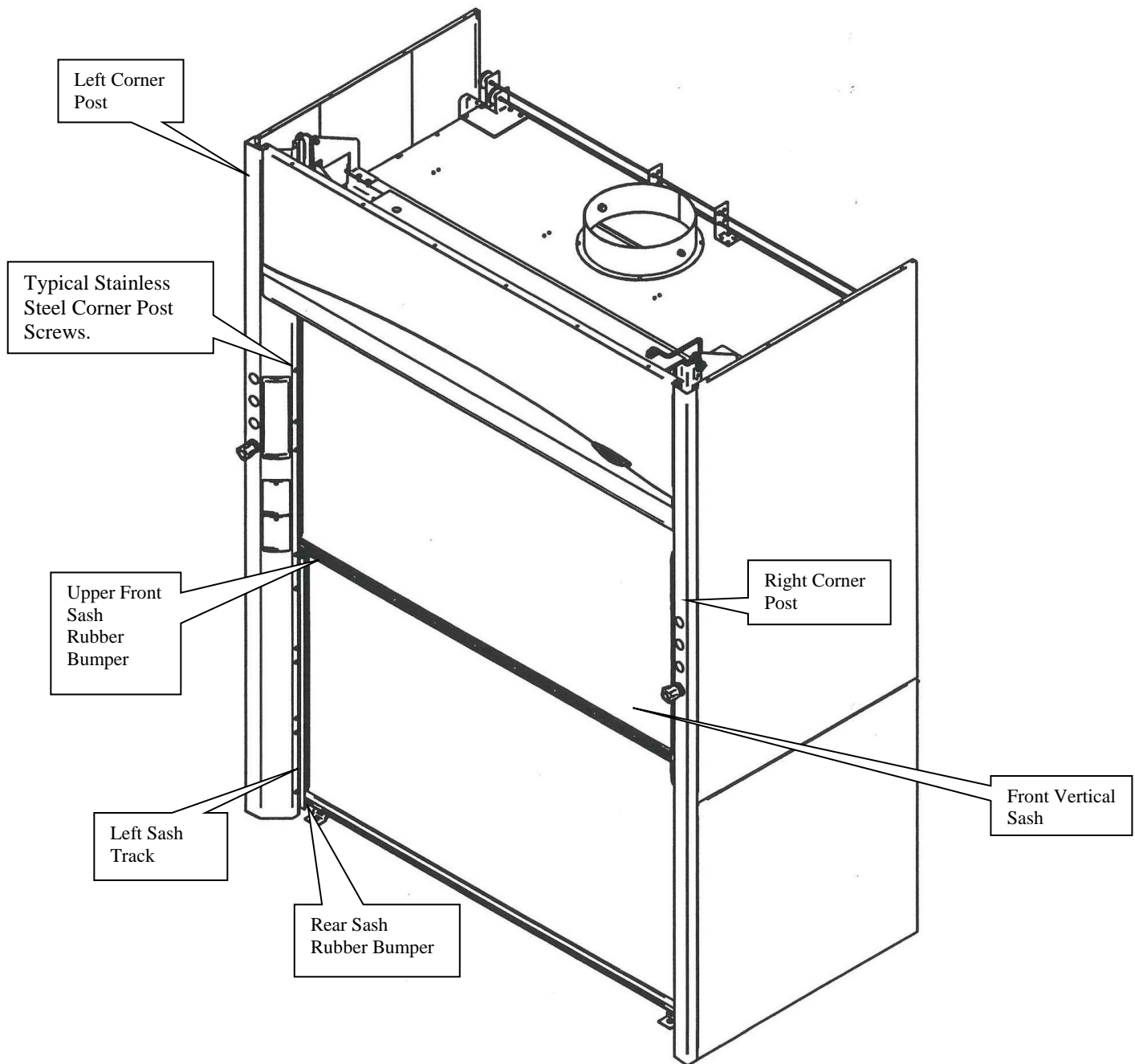
Figure 3-3

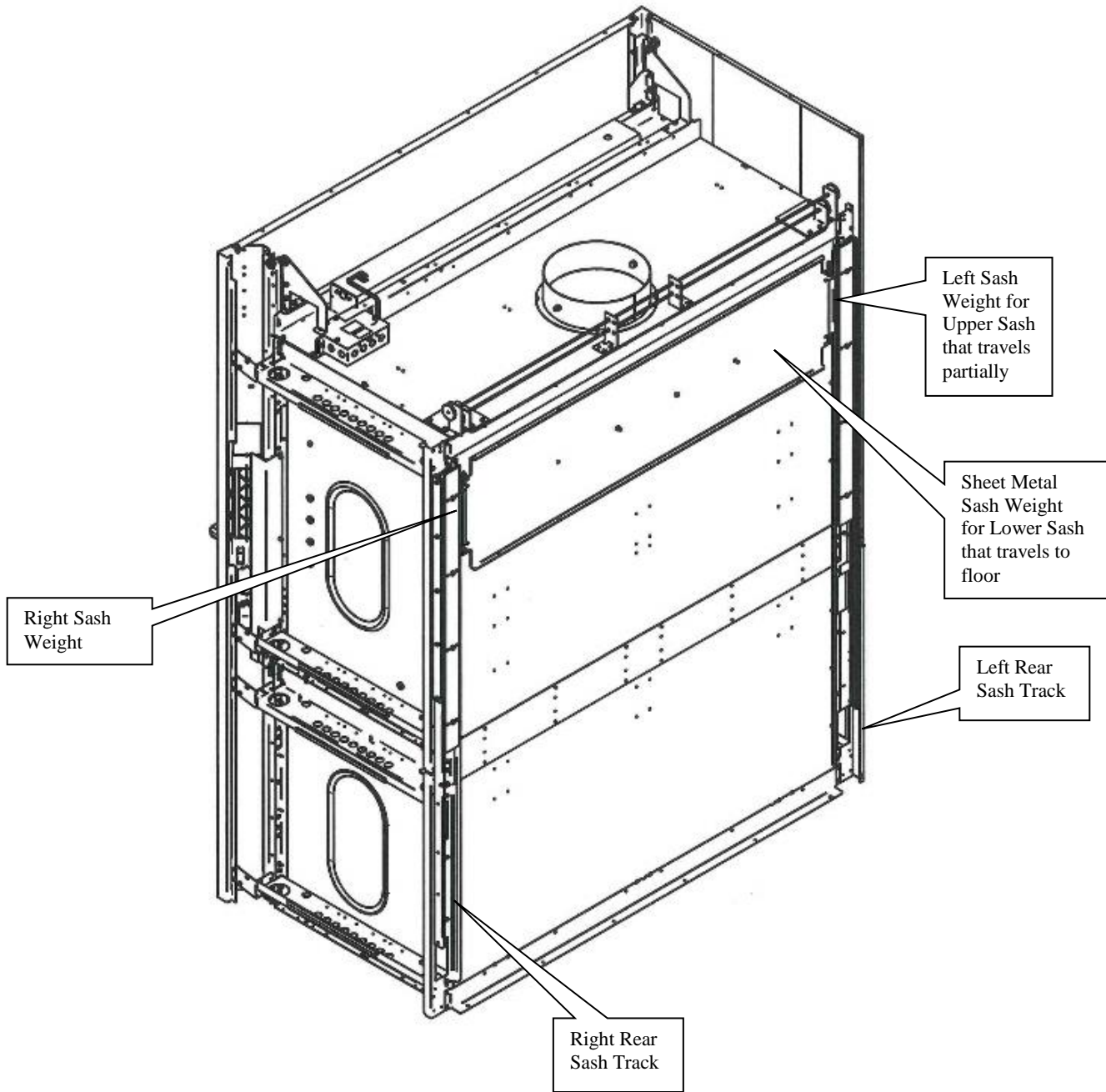Rear View

Figure 3-4

Lower Sash Track & Threshold Installation for Horizontal-Sliding Sashes on 8', 10', 12', & 16' Floor-Mounted Hoods

H

The lower sash track assembly is comprised of a sash track threshold, front and rear ramps, and threshold support brackets. With the corner posts removed the sash track threshold should be placed on the floor between the left and right corner posts, with their corner posts removed at this time. Lay the threshold flat, perpendicular to the hood liner walls. The sash track threshold will extend to the hood sidetracks. Once the sash track threshold has been placed in its final position, check to be sure it is level. It must be level to ensure proper movement of each of the door assemblies. If the threshold is not level, place shim material under it until it becomes level. Attach the sash track threshold to the hood sidetracks with the threshold support brackets and #10-24 hardware. Both front and rear ramps are held in position without mechanical fasteners and can be removed for easy cleaning, should it be required. Simply place the flange of each ramp assembly in the position as shown in the diagram below for proper fit. The ramps are self-leveling to match your existing floor conditions. The hood corner posts should be installed prior to the placement of the front sash ramp. See Figure 3-5 and Figure 1-2.


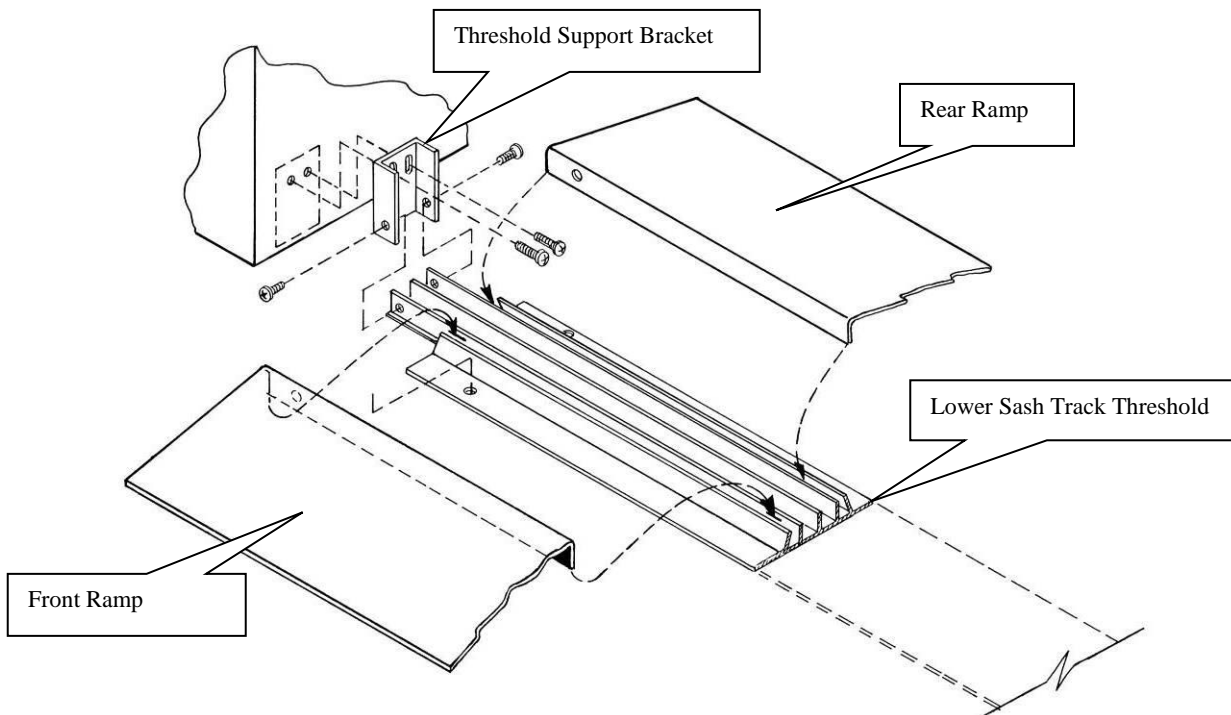
Figure 3-5

H Horizontal-Sliding Sash Installation for the 8', 10', 12', & 16' Floor-Mounted Hoods

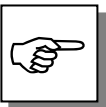The Floor-Mounted Hood with horizontal-sliding sashes is supplied with four individual sash doors. These doors have been packaged separately during shipment and require installation in the field. The doors should not be installed until the lower sash track threshold has been properly installed. Each door assembly is provided with two mounting brackets with nylon roller bearings. The roller bearings have been positioned in the center location of each mounting bracket to provide for maximum door adjustment. The upper header sash track features two tracks that will support two of the door assemblies. Raise the door panel up and into the upper track section on your hood and place the nylon roller bearings directly in the track section. At the same time, position the bottom edge of the door assembly into the corresponding lower track slot. As each door is placed into position, slide the door the full length of the track assembly to assure proper height adjustment. If the door drags in the lower track, adjust the nylon roller bearing on the mounting bracket to raise the height of the door. Reverse the movement of the bearing should the door assembly be too short and not fit in the lower track properly. An average space of 1/4" to 5/16" should be left between the bottom of each door and the bottom of the lower sash track for proper door travel. Repeat this procedure for the four door assemblies and place two of the doors on each of the front and rear track runners. Once proper door operation has been achieved, the sash track threshold should be permanently attached to the floor. Anchor bolts should be used in the holes provided in the lower sash track threshold mounting. The anchor bolts are not supplied. RTV sealant can also be used to hold the lower sash track threshold in position during normal operation. See Figure 3-6 and Figure 1-2.

TIP: SHOULD YOU EXPERIENCE SLIGHT BINDING OF THE SASH DOORS OR DOOR IN THE TRACK, SPRAY THE TWO SASH GLIDES ON THE BOTTOM OF THE DOOR FRAME WITH SILICONE SPRAY.

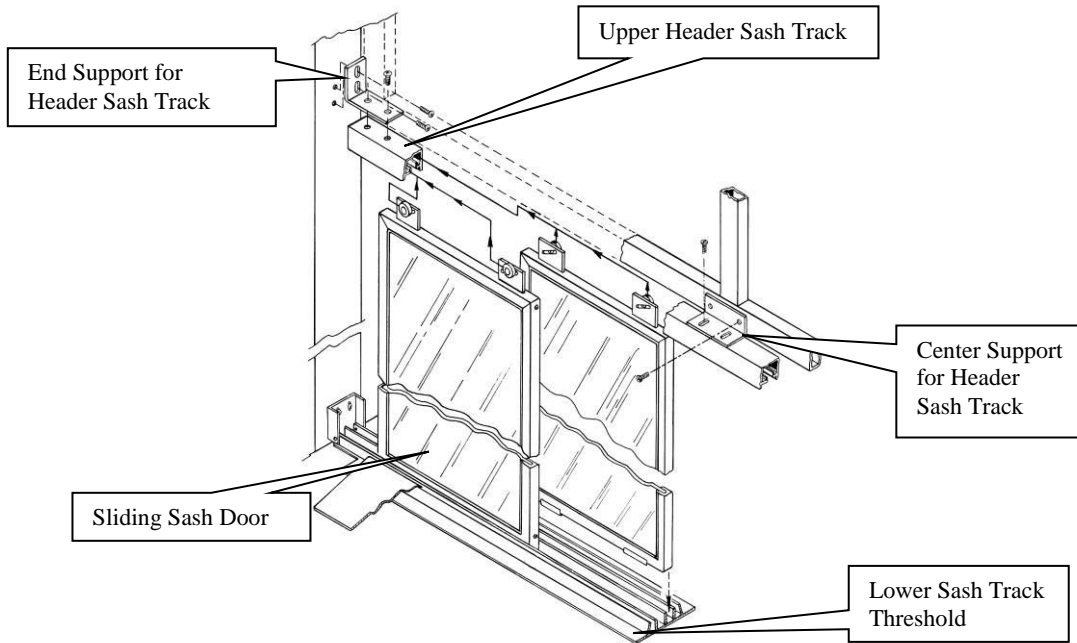
Figure 3-6

Corner Post Installation for the Floor-Mounted Hoods

Both left and right corner posts have been shipped uninstalled. The one-piece corner posts are to be installed once the upper and lower cabinet assemblies have been properly positioned.

The edges on the corner posts fit directly onto the side frames. The front inner edge of both corner posts are held in place by stainless steel machine screws. The outer back edge of both corner posts are held in place with steel self-drilling screws. The screws are included in the hood manual packet. See Figure 3-2 and Figure 3-3.

Baffle Installation for the Floor-Mounted Hoods

The baffles are critical to the proper airflow and performance of the hood. Be sure the baffles are resting in the proper baffle mount supports. See Figure 3-7 for various sizes of baffles and their orientation.

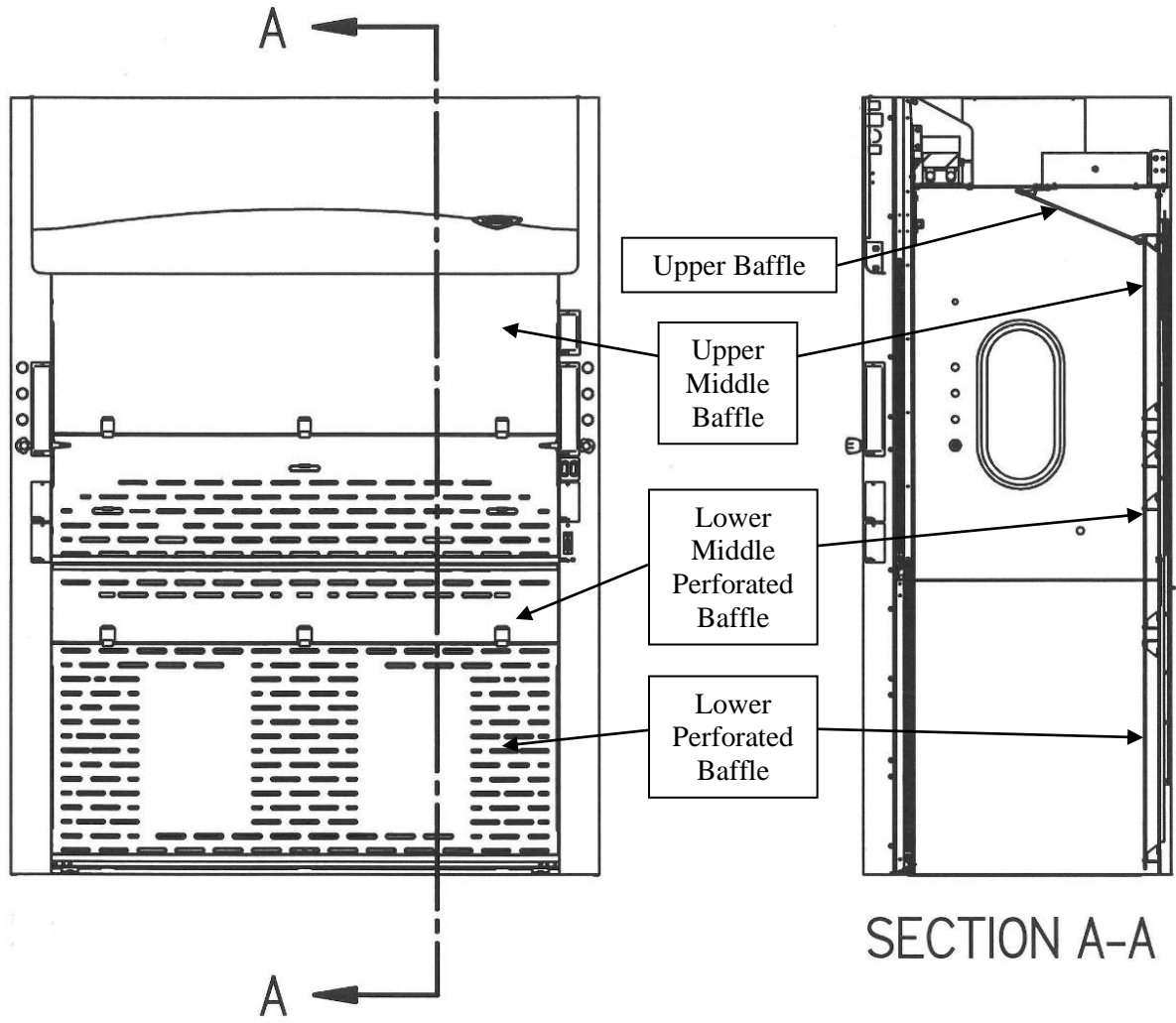
Figure 3-7

Header Installation for the Floor-Mounted Hoods

The header is shipped separately in its protective packaging. To install, the header is fastened to the corner covers by four #12 screws. Reach behind the corner posts from the side and install the screws to support the header. See Figure 3-8.

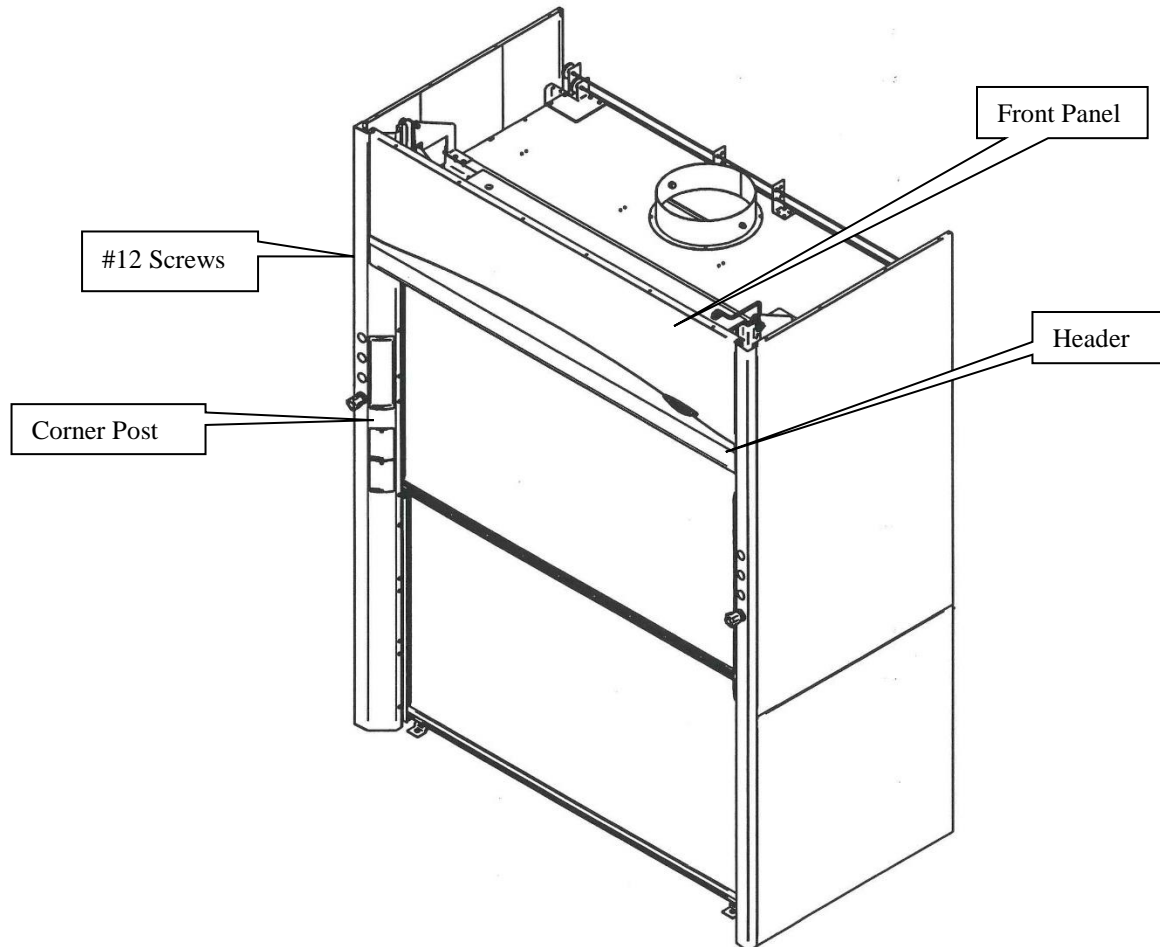
Figure 3-8

Front Panel Installation for the Floor-Mounted Hoods

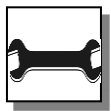
The hood front panel(s) is shipped separately and is protected. To install each panel, hang the two plastic cylinders on the backside of the front panel over the corner posts and center support (10', 12', 16' models). The bottom of the front panel will then slip behind the header once it has been properly secured at the top. See Figure 3-3 and Figure 3-8.

Connecting to the Hood Exhaust System

WARNING: The weight of the exhaust ductwork system must be supported independently of the hood superstructure. Do not allow this weight to be supported by the hood structure as damage to the hood may occur.

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The exhaust connection should be installed by a qualified HVAC contractor. The exhaust connection on your hood has been designed for 12" nominal pipe (12.75" OD) to allow for minimum static pressure loss with proper transport velocities away from the hood. Consult Labconco Customer Service should you require help sizing your blower for the exhaust volume and total system static pressure loss.

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The selected exhaust duct material should match the hood procedures and chemicals used to ensure compatibility.

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Connecting the Electrical Supply Source to the Protector Fume Hood

Prior to connecting any electrical wiring to the fume hood structure, refer to the hood identification plate for the proper electrical requirements of your specific model.

WARNING: The building electrical supply system for Protector Hoods should include overload protection. A switch or circuit breaker should be in close proximity to the equipment and within easy reach of the operator. The switch or circuit breaker is to be marked as the disconnecting device for the equipment. Consult the NEC-2002 for proper installation.

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The identification plate, model number, serial number, and electrical connection boxes are accessible from the front of the fume hood by removing the front panel.

The Protector XL Hood is normally wired for 115 Volt, 50/60 Hz, 20 Amp or 230Volt, 50/60 Hz, 10 Amp electrical service. Check the I.D. plate behind the front panel for voltage verification. The number of circuits varies depending on the model. All of the electrical connections are terminated at the field wiring terminal box for hook-up by a qualified electrician. We recommend each circuit be a dedicated branch circuit. However, if wired together the maximum load allowed is the sum of individual outlets plus the rating of the unit (i.e. 2 Amps). The single point internal junction box is used for the connection of the lights, blower, and duplex outlets. Refer to the wiring diagram for your Protector XL in *Appendix C: Protector XL Specifications*.

The fume hood is required to be grounded to the MAINS protective earthing ground for safe operation. Using a ring terminal sized for a 10-24 machine screw, connect the MAINS ground conductor to the grounding lug marked with the protective earthing symbol, . Only MAINS ground conductors should be connected to the protective earthing ground lug, no other conductors should be connected to this grounding lug. Using wire nuts connect the MAINS supply conductors to the fume hood supply wires. Insure that the wires are connected as per the appropriate wire color codes for the input voltage. For 115V Phase (Hot) is black and Neutral is white, for 230V Phase1 is brown and Phase2 is blue. Refer to the wiring diagram for your Protector XL in *Appendix C: Protector XL Specifications*.


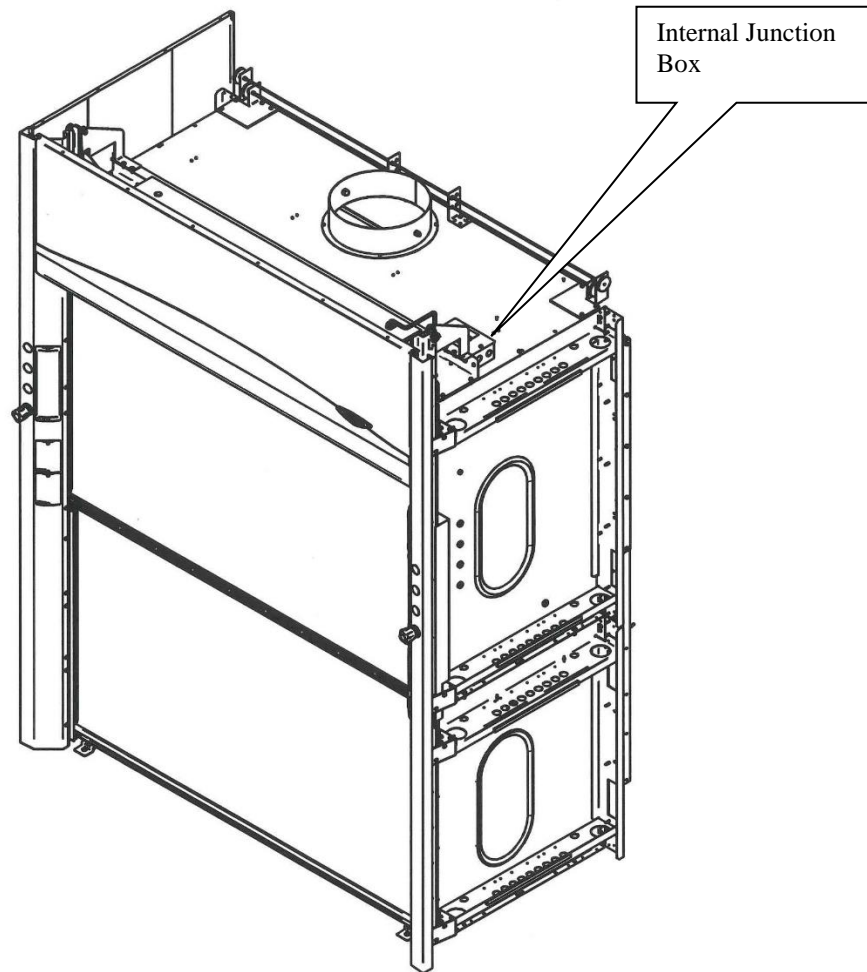
Figure 3-9

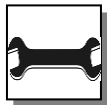All wiring for the fume hood **SHOULD** be performed by a licensed electrician and conform to all local codes. In most cases, the hood will require the use of shielded conduit to protect the wiring into the hood. The grounding connection shall not be made to the terminal box cover.

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The LED light has been mounted outside the top liner panel and is sealed from vapors inside the hood structure. To change the fluorescent light bulbs in your hood, you must first remove the front panel from the hood. Next remove the knock out plugs holding the light fixture in place. Lift fixture up and replace any defective bulbs. Reverse order to reassemble.

Connecting the Service Lines to the Protector Fume Hood

The hoods with service fixtures have been plumbed from the valve to the hose connector or gooseneck for your installation convenience. Supply tubing shall be provided by the qualified installer. Tubing can enter the hood from above, through the back, or through the work surface to make these connections to the service fixtures.

NOTE: Inspect all fittings for leakage. Tighten the fittings slightly if needed.

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CAUTION: Do not use oxygen with any standard service fixture. Contact Labconco Customer Service for oxygen fixture information.

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Should access to the hood plumbing fixture bodies be required, remove the service access plate on the hood front corner posts by loosening their individual screws (Figure A-1 in *Appendix A*). The valve body will now be fully exposed for any service work that may be necessary. The service fixtures supplied on your laboratory hood are designed for use with the following services:

- Air
- Hot Water
- Vacuum
- Cold Water
- Natural Gas – See caution below

WARNING: Contact Labconco Customer Service directly before using any service other than those listed above in these valves to assure full compatibility.

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CAUTION: Natural gas should be used only in the service fixture that has been pre-plumbed with brass tubing. Sulfur content of the gas could cause deterioration of standard copper supply lines.

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Sealing the Protector Hood to the Floor

When the hood has been set in place, ducted, wired, and plumbed, it should be sealed at the floor to prevent spilled materials from collecting under the walls of the hood. Materials such as silicone sealants are recommended to seal the hood structure to the floor.

Certifying the Protector Fume Hood

The combination of your laboratory hood, exhaust ductwork, and exhaust blower gives you the flexibility to change the airflow at the sash opening of your hood. To determine the actual face velocity at the sash opening, airflow velocity readings will need to be taken. This should be done across the sash opening of the hood in accordance with the *Industrial Ventilation Manual* section on laboratory hoods (see *Appendix E: References*). Labconco recommends an average face velocity at the sash opening of 60 to 100 feet per minute. Consult Chapter 2 for proper airflow volumes for your particular model.

Your Protector Fume Hood has been tested at the factory per ASHRAE 110-1995. All hoods achieve an “as manufactured rating” of less than 0.05 part per million (ppm) at 4 liters per minute (lpm); AM<0.05 (consult Labconco for individual fume hood ratings). For “field use” ASHRAE testing contact Labconco Sales Engineering Team or Customer Service for a certified on-site contractor.

NOTE: Face velocity profiles and smoke testing should be done periodically to ensure safe performance.

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CHAPTER 4

PERFORMANCE FEATURES AND SAFETY PRECAUTIONS

Performance Features:

The Protector XL Floor-Mounted Laboratory Hood is designed to meet the needs of the laboratory scientist who must transport equipment and products for experimentation and testing at OSHA approved high performance/low flow velocities for vertical sashes as low as 60 feet per minute and horizontal sashes as low as 80 feet per minutes. Vertical Floor-Mount Hoods are high performance. The laboratory fume hood has been designed to effectively contain toxic, noxious, or other harmful materials when properly installed. A fully featured by-pass hood with perforated baffle and by-pass block, this hood maintains safe airflow. Optional energy saving A-Style Combination Sash models are also available. The hood features by-pass airflow design that promotes full containment as the sash is moved. Airflow is diverted behind the front panel and through the perforated sash handle to help control fluctuations in face velocity, which occur as the sash is closed.

- a. Vertical:** Unique sash provides maximum visibility of 73.5" high while conserving energy by limiting sash travel to 28". Vertical-rising sash may be raised from a closed to 64" operating height. The fume hood features a containment-enhancing sash handle. Vertical Floor-Mount Hoods are high performance. Optional sash stops are available to limit sash height and reduce energy usage.
 - b. Horizontal:** The horizontal sliding sash doors provide maximum visibility and load height of 73.5". The total exhaust volume is based on half the opening on the 8' model and one quarter the opening on 10', 12', and 16'.
- By-pass airflow design** ensures containment as the sash moves.

Chapter 4: Performance Features & Safety Precautions

3. **Large usable interior work depth and interior height of 84"** provides ample working space.
4. **Baffle** directs airflow to the rear of the interior to provide efficient airflow. The baffle may be removed for cleaning purposes only. The Opti-Zone™ baffle provides optimized airflow and face velocity profile for optimal containment.
5. **Exterior access cover plates** are removable for easy access to plumbing valves and sash adjustment hardware when access through the sides is not available.
6. **Lift-Away™ front panel** provides easy access to electrical wiring, sash weights, and lighting fixtures.
7. **Energy efficient LED lighting** is located behind a laminated safety glass shield mounted to the top of the hood. The factory-wired instant start T8 lighting is serviceable from outside the hood cavity. Additionally, the long lasting 50,000 hour direct wired LED T8 bulbs are more energy efficient, utilizing approximately ½ the power of fluorescent bulbs. See Appendix A for wattage.
8. **Low mounted, factory-wired light and blower switches** are ADA compliant.
9. **Streamlined corner posts** provide maximum visibility and the flexibility to add services after installation.
10. **All hoods are factory prepared for up to 8 service fixtures.**
11. **Duplex electrical receptacles** are mounted on the right and left corner posts as requested. Receptacles are factory-wired to hood single point junction box.
12. **Accessory Guardian™ Digital Airflow Monitor or Guardian Airflow Monitor** continuously monitors face velocity. An audio/visual alarm alerts the user to low airflow conditions. The right corner post is factory prepared to accommodate the Guardian Monitor (sold separately).
13. **Optional Sash Models.** 4'-8' Vertical Sashes are offered in smooth anti-racking cable. 8'-16' Horizontal Sashes slide laterally to open. For other sash options, contract Labconco.
14. **Frame of epoxy-coated steel and aluminum** is durable and corrosion resistant.
15. **Exhaust Connection.** The hood features 12" (12.75" OD pipe) exhaust connections sized to allow for a minimum static pressure loss through the hood structure while providing a good transport velocity through the exhaust system.
16. **Optional Ceiling Enclosure Kits** are available for a decorative facade between the hood and the ceiling.
17. **Optional Sash Stops** provide a means of controlling the operating height of the upper vertical sash.

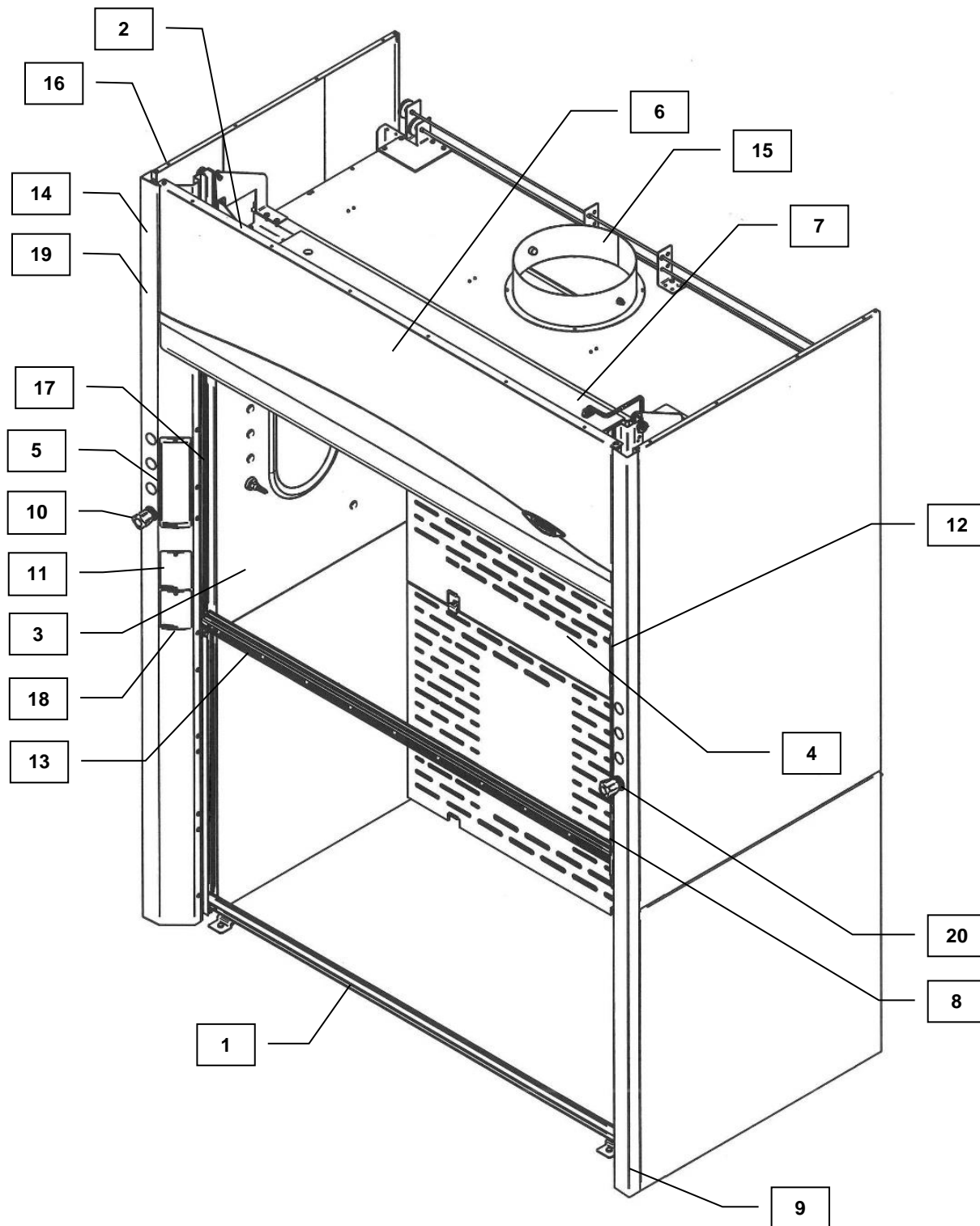
Figure 4-1

Chapter 4: Performance Features & Safety Precautions

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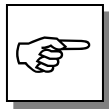
Safety Precautions

Although the laboratory hood has been engineered to maintain optimum operator safety, caution should always be used while working in the hood. Prior to using the hood, check to make sure that the exhaust blower is operating and that air is entering the hood at its specified face velocity.

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USE GOOD HOUSEKEEPING IN THE HOOD AT ALL TIMES. CLEAN UP SPILLS IMMEDIATELY WITH A MILD DETERGENT. PERIODICALLY CLEAN HOOD INTERIOR, INCLUDING FLUORESCENT LIGHT GLASS PANEL. REPLACE BURNED OUT LIGHT BULBS TO MAINTAIN MAXIMUM ILLUMINATION.

DO NOT OVERLOAD THE WORK SURFACE WITH APPARATUS OR WORK MATERIAL. THE SAFE OPERATION OF THE LABORATORY HOOD IS BASED UPON HAVING PROPER AIRFLOW THROUGH THE STRUCTURE. DO NOT PLACE LARGE, BULKY OBJECTS SUCH AS BLOCK HEATERS, DIRECTLY ON THE HOOD WORK SURFACE. INSTEAD, ELEVATE THE OBJECT 2" TO 3" ON BLOCKS TO ALLOW A FLOW OF AIR UNDER THE OBJECT AND INTO THE LOWER REAR BAFFLE EXHAUST SLOT. ENSURE BLOCKS ARE LEVEL AND SECURED IN PLACE.

Blocking the bottom of the baffle at rear of hood will change the airflow pattern in the hood causing turbulence and possible leakage at the face of the hood. (Don't store containers or supplies against baffles, as this will affect airflow through the hood).

Avoid placing your head inside hood. Keep hands out of hood as much as possible.

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For Vertical Sashes, always work with the lower sash closed. For Horizontal Sash Door models, always work with a maximum of two doors open for 8' models and one door open for 10', 12', and 16' models. Refer to Chapter 2 for airflows.

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Always work as far back in hood as possible. It is best to keep all chemicals and apparatus 6" inside the front of the hood.

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This hood does not feature explosion-proof electrical components, unless ordered separately. Therefore, use of flammable or explosive materials in quantities above the explosive limit are not recommended.

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Do not work with chemicals in this hood without the exhaust system running. Do not store chemicals in a fume hood.

Chapter 4: Performance Features & Safety Precautions

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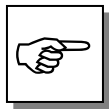
Perchloric acid use in this hood is prohibited.

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High level radioisotope materials are prohibited for usage in this hood.

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AVOID CROSS DRAFTS AND LIMIT TRAFFIC IN FRONT OF THE HOOD. AIR DISTURBANCES CREATED MAY DRAW FUMES OUT OF THE HOOD.

The use of heat-generating equipment in this hood without the exhaust system operating properly can cause damage to the hood.

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The Protector Laboratory Hood should be certified by a qualified certification technician before it is initially used. The hood should be re-certified whenever it is relocated, serviced or at least annually thereafter.

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Ensure that the unit is connected to electrical service in accordance with local and national electrical codes. Failure to do so may create a fire or electrical hazard. Do not remove or service any electrical components without first disconnecting the hood from electrical service.

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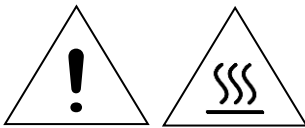
Proper operation of the fume hood depends largely upon the hood's location and the operator's work habits. Consult the Reference Manual in *Appendix E*.

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If the unit is not operated as specified in this manual it may impair the protection provided by the unit.

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Do not touch the blower motor. The surfaces of the motor can become hot and could cause burns.

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Do not position the fume hood so that it is difficult to operate the main disconnect device.

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To prevent the possibility of minor injury keep hands and fingers clear of sprockets at the top of the four corners.

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CHAPTER 5

USING YOUR PROTECTOR XL

S Operating the Vertical-Rising Sash

4' **5'** Because of the Protector Hood counterbalanced sash mechanism, it will take only a few pounds of force to move the sash up or down, and you can operate the sash smoothly with one or two hands positioned anywhere along the handle. The vertical-rising sash may be raised to a maximum 64" operating height. The airflow requirements should be sized for the 64" operating height; if using sash stops then the airflow requirements can be reduced by approximately 40% at 54". See *Chapter 2* for airflows.

6' **8'**

H Operating the Horizontal-Sliding Sashes

8' **10'** Horizontal-sliding sashes allow the operator the flexibility of arranging the sashes at a position suitable for various procedures. The glass sashes have a 73.5" sash opening providing a large viewing area. The airflow requirements are sized for the 50% open sash condition for the 8' model and 25% for the 10', 12', and 16' models.

12' **16'**

A Operating the A-Style Combination Sash

Some other hood models have additional energy saving sashes called A-Style Combination Sashes in place of vertical-rising sashes. These combination sashes allow the operator to use the hood with sashes either half open horizontally or vertically to conserve energy. The horizontal sashes are used in normal operating mode. Optional sets of sash stops can be installed to prevent raising the vertical sash above the half-open and fully-closed positions unless manually defeated by the operator. The airflow requirements are sized for the 50% open sash condition.

S **A** **Operating the Blower**

Your Protector Fume Hood utilizes a remote style blower, which can be activated by turning the blower switch to “ON.” You can validate the hood performance by watching smoke drawn into the hood face opening.

S **A** **Operating the Lights**

Your Protector Fume Hood utilizes a factory-wired LED light to illuminate the hood interior (see Appendix A). Simply turn the light switch to “ON” to operate.

Working in your Protector Fume Hood

Planning

- Thoroughly understand procedures and equipment required before beginning work.
- Arrange for minimal disruptions, such as room traffic or entry into the room while the hood is in use.

Start-up

Labconco follows the recommendations of the American National Standard Institute with regards to fume hood exhaust systems. A full list of requirements can be found in the ANSI Standard No. Z9.5 *Laboratory Ventilation*. If you are unsure of the proper use of the exhaust blower for your fume hood, contact your safety officer or industrial hygienist before making any modifications to the blower’s operational status (on/off/low, etc.). Keeping the exhaust system under constant negative pressure is an excellent way to ensure hazardous fumes in the fume hood do not enter the laboratory space inadvertently.

- Turn on the light.
- Slowly raise the sash and verify the fume hood has proper airflow.
- Check the baffle air slots for obstructions.
- Allow the hood to operate unobstructed for two minutes.
- Wear a long sleeved lab coat and rubber gloves. Use protective eyewear. Wear a protective mask if appropriate.

Loading Materials and Equipment

- Only load the materials required for the procedure. Do not overload the hood.
- Do not obstruct the front air foil (sill), or rear baffle slots.
- Large objects should not be placed close together and spaced above the work surface to permit airflow to sweep under the equipment.
- After loading the hood, wait one minute to purge airborne contaminants from the work area.

Chapter 5: Using Your Protector XL

Work Techniques

- Keep all materials at least 6 inches inside of the sash and at least two inches from the side walls and perform all contaminated operations as far to the rear of the work area as possible.
- Keep all clean and contaminated materials in the work area separate for clear identification.
- Avoid using techniques or procedures that disrupt the airflow patterns of the hood.

Final Purging

- Upon completion of the work, the hood should be allowed to operate for two minutes undisturbed, to purge airborne contaminants from the work area before closing the sash.

Unloading Materials and Equipment

- Objects in contact with contaminated material should be surface decontaminated before removal from the hood.
- All open trays or containers should be covered before being removed from the hood.

Shutdown

- Turn off the light and close the sash.

CHAPTER 6

MAINTAINING YOUR PROTECTOR XL

Now that you have an understanding of how to work in the fume hood, we will review the suggested maintenance schedule and the common service operations necessary to maintain your fume hood for peak performance.

Service Safety Precautions

- If performing any electrical maintenance, always disconnect the power at the main disconnect.
- If performing decontamination inside the fume hood, consult your safety officer for proper personal protective equipment and procedure.
- Since some service operations require a step ladder, always use proper safety and consult your safety officer
- If performing maintenance on any service lines, always shut off the supply first.
- Some removable components may be heavy, follow safe-lifting guidelines.
- Verify all components are installed correctly with performance verified before conducting normal operations.

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Chapter 6: Maintaining Your Protector XL

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Routine Maintenance Schedule

Weekly

- Using ordinary dish soap to clean the surface inside of the fume hood.
- Using an appropriate glass cleaner, clean the sash and all glass surfaces.
- Operate the fume hood blower, noting the airflow velocity through the hood using a source of visible smoke.

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Biannually

- The sash assembly should be checked to ensure proper operation and to make sure there are no signs of abnormal wear on the sash pulleys, cables and clamps.

Routine Service Operations

Front Panel Removal:

1. Simply lift the front panel up and then away from the hood to provide access to the top.


Changing the Lamp:

1. Turn light switch to “OFF”.
2. Remove the front panel as noted earlier.
3. Reach over the front header of the hood and remove knock out plugs at both ends of fixture. Lift fixture up.
4. Remove the lamp by pushing it out of the spring-loaded lamp socket and swinging it out of the other lamp socket.
5. Install the new lamp by reversing the removal procedure.

CHAPTER 7

MODIFYING YOUR PROTECTOR XL

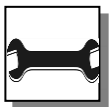There are several ways to modify the fume hood for your individual requirements. These include the addition of work surfaces, service fixtures, air monitor, distillation grids, electrical duplex outlets, and ceiling enclosures.

Installing Work Surfaces

Contact Labconco Customer Service for ordering information of optional work surfaces.

Installing Ceiling Enclosures Above the Fume Hood

Your Protector Fume Hood has mounting holes to accept a ceiling enclosure to close off the area between the top of the hood and the ceiling. Contact Labconco Customer Service for ordering information. Labconco offers both non-adjustable fixed height ceiling enclosures as well as adjustable ceiling enclosures per the charts below. Other special sizes are available upon request.

3-Sided Ceiling Enclosure Kits

The panels extend above the top of the hood to the ceiling to hide exposed ductwork, plumbing and wiring. For any 3 ft to 8 ft Protector XL Benchtop, Premier, XStream, Stainless Steel or PVC Fume Hood. Each kit includes three glacier white powder-coated steel ceiling enclosure panels, one for the front of the hood and one for each side of the hood.

Configure Your Non-Adjustable Fixed Height Ceiling Enclosure Kit Catalog Number

Use this key to configure the seven digit catalog number to order your 3-Sided **Fixed Height** Ceiling Enclosure Kit. For example, **955011** is a 3-Sided Ceiling Enclosure Kit for a 5' x 37.7" deep Protector XL Floor-Mounted Hood with a fixed height of 12".

9 5 5 0

STEP 1. Select the **nominal width** of your fume hood. This number is the fifth digit of your catalog number.
 4 = 4'
 5 = 5'
 6 = 6'
 8 = 8'

STEP 2. Select the **depth** based on the model chosen. Choose 0 for Premier and 1 for XStream. Choose 1 for PVC. Choose 0, 1 or 2 for the correct depth of the XL Benchtop model. Choose 0 or 1 for the correct depth of Stainless Steel. This is the sixth digit of your catalog number.

No.	Ext. Depth	Model
1	37.7"	XL Floor-Mounted
2	43.7"	XL Floor-Mounted

STEP 3. Choose the correct **height** that matches the number. This height should match the height between the top of the hood and the building ceiling. This is the seventh digit of your catalog number.

No.	Height
0	6"
1	12"
2	18"
3	24"
4	30"
5	36"

Chapter 7: Modifying Your Protector XL

Configure Your Adjustable Height Ceiling Enclosure Kit Catalog Number

Use this key to configure the seven digit catalog number to order your 3-Sided **Adjustable Height** Ceiling Enclosure Kit. For example, **9425511** is a 3-Sided Ceiling Enclosure Kit for a 5' x 37.7" deep Protector XL Floor-Mounted Hood with a height range of 14-19".

9**4****2****5**

STEP 1. Select the **nominal width** of your fume hood. This number is the fifth digit of your catalog number.

4 = 4'

5 = 5'

6 = 6'

8 = 8'

STEP 2. Select the **depth** based on the model chosen. Choose 0 for Premier and 1 for XStream. Choose 1 for PVC. Choose 0, 1 or 2 for the correct depth of the XL Floor-Mounted model. Choose 0 or 1 for the correct depth of Stainless Steel. This is the sixth digit of your catalog number. *Order 9548700 separately.

No.	Ext. Depth	Model
1	37.7"	XL Floor-Mounted
2	43.7"	XL Floor-Mounted

STEP 3. Choose the correct **height range** that matches the number. This height range should match the height between the top of the hood and the building ceiling. This is the seventh digit of your catalog number.

No.	Height Range
0	11-14"
1	14-18.6"
2	18.6-24.4"

Note: The previous generation listed as 9414-W-D-H is presented below to aid product service.

9**4****1****4****W****D****H**

Note: *For previous generation Protector XL Floor-Mounted Hoods, Ceiling Enclosure Conversion Kit #9548700 is required for unique rear mounting brackets and must be ordered separately.

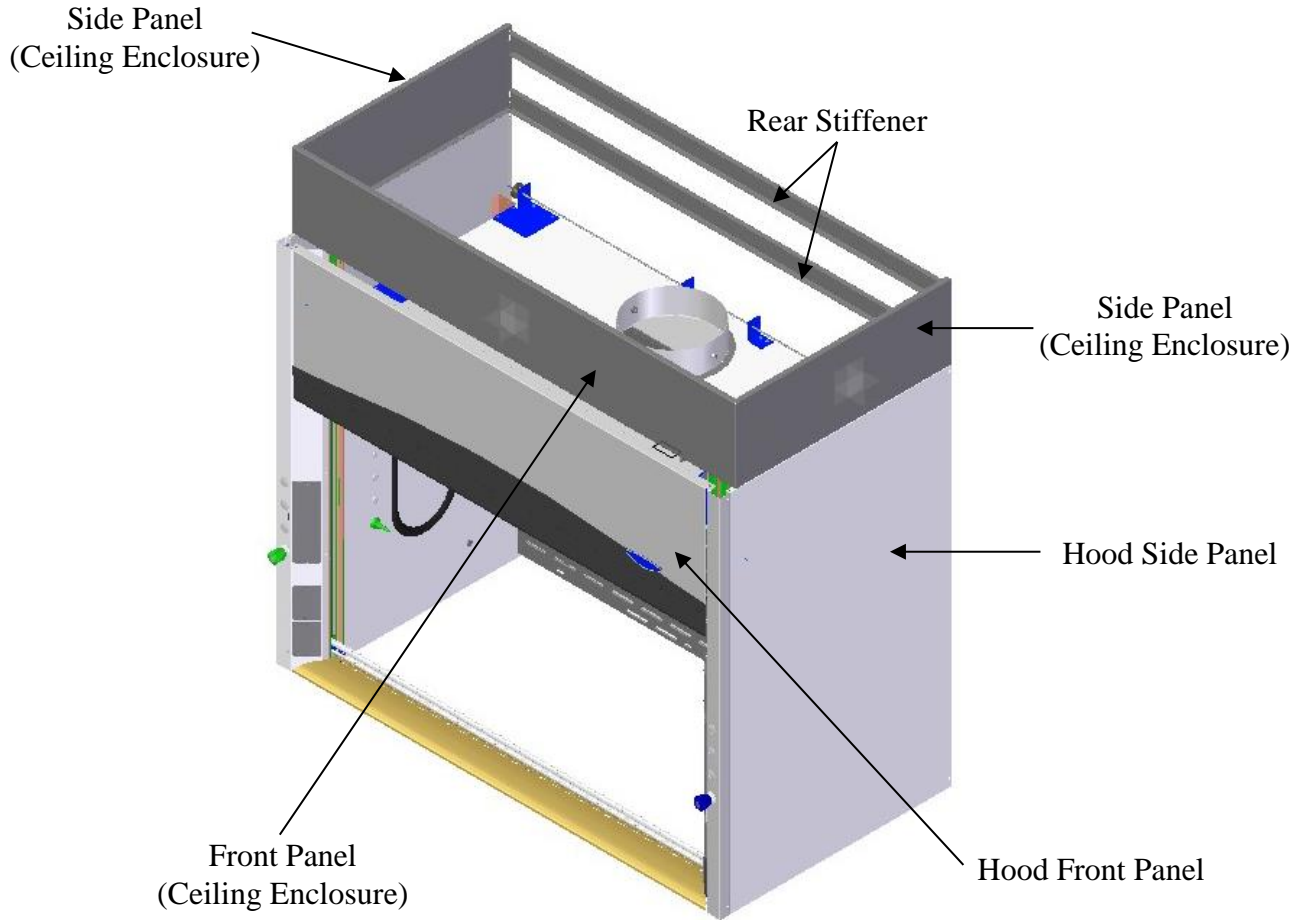
Figure 7-1

Installing Additional Service Fixtures

Additional service fixtures can be installed in the available service fixture holes in both sidewalls and corner posts. The fume hood is factory set to accept up to four valves per side. Contact Labconco Customer Service for information.

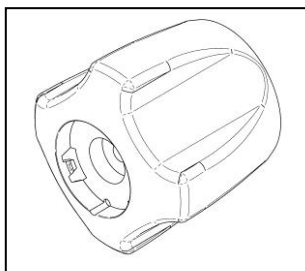
Figure 7-2 - Knob

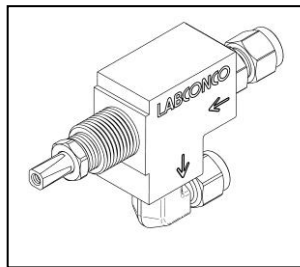
Figure 7-3 - Valve

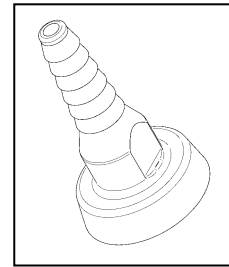
Figure 7-4 - Hose Connector

Chapter 7: Modifying Your Protector XL

Installing Guardian™ Digital Airflow Monitor or Guardian™ Airflow Monitor

The Guardian Airflow Monitors continuously monitor face velocity through the fume hood opening. The fume hood right corner post is factory prepared to mount either monitor. Contact Labconco Customer Service to order.

Floor Mount Distillation Grids – Field Installation

The floor mount distillation grids have been strategically placed with the vertical rod center lines in front of the lower baffle and middle baffle. The floor mount distillation grids allow the hood user to mount glassware, motors, stirrers, and other apparatus. Contact Labconco Customer Service for ordering information for Floor-Mounted Hoods.

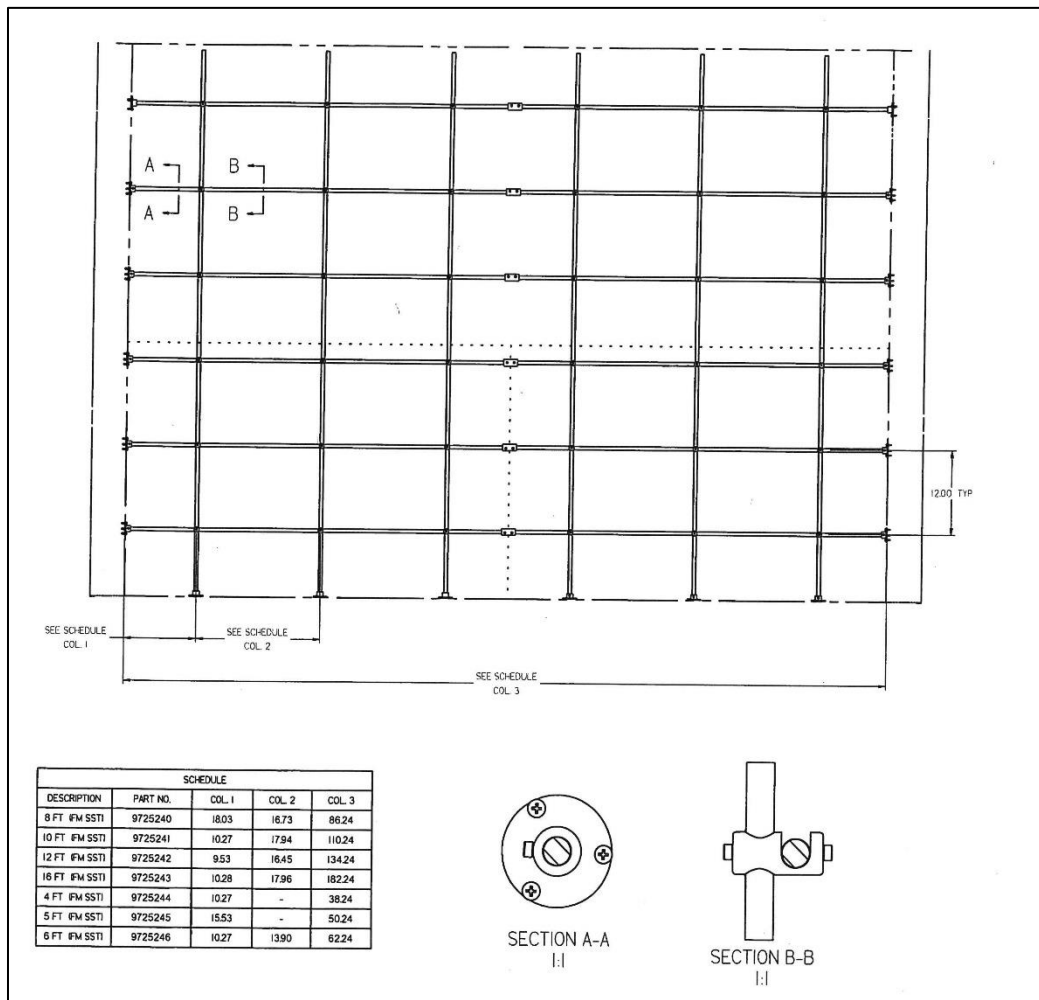
Figure 7-5

Sash Stop Kit – Field Installation P/N 9410300

The sash stop kit restricts how far a vertical-rising sash may be opened. This small plastic device may be easily field installed on the fixture corner post of any fume hood.

Installing an Electrical Duplex Outlet

Your Protector Fume Hood can be ordered with duplex outlets, however, if you ordered a model without an electrical duplex outlet you can have one installed in the field by a qualified electrician. Contact Labconco Customer Service for ordering information. (Not acceptable on explosion-proof hoods).

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
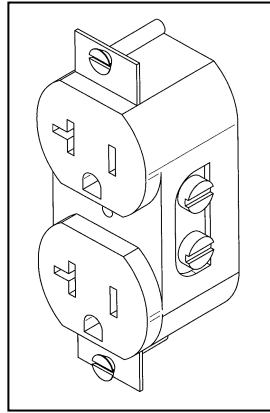
Figure 7-6

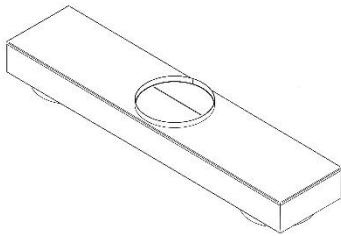8' Dual Exhaust Adapters – Field Installation

The Type 304 stainless steel single exhaust adapter allows 8' Protector Hoods to be exhausted from two fume hood exhaust collars spaced 48" apart to a single 12" (12.81 ID) exhaust, centered on the hood. The single centered exhaust outlet is located 9.3" from the rear of the hood for all 8' Protector Hoods.

For other special sizes, consult Labconco. The performance specifications for the 8' Dual Exhaust Adapters are listed below.

Catalog No.	Description	Dimensions (w x d x h)	Height Above Hood (inches)	Equivalent Resistance (feet)	Shipping Wt. (lbs.)
9595108	Dual Exhaust 12 x 12 x 12 x 6 High Flush	63" x 14.2" x 8.5"	0	150	60
9595109	Dual Exhaust 12 x 12 x 12 x 9 High	63" x 14.2" x 11.5"	3	120	70

CHAPTER 8

TROUBLESHOOTING

Refer to the following table if your fume hood fails to operate properly. If the suggested corrective actions do not solve your problem, contact Labconco for additional assistance.

PROBLEM	CAUSE	CORRECTIVE ACTION
Remote blower and lights won't operate	Wires not connected at junction boxes or switches.	Check connection of switches. Check connection to control box on top of unit.
	Circuit breakers tripped in building electrical supply.	Reset circuit breakers.
Remote blower won't operate, but lights work	Blower wiring is disconnected.	Inspect blower wiring and switch.
	Belt broken	Replace belt.
	Blower motor is defective.	Replace blower motor.
Fume hood blower operates but lights will not operate	Lamp not installed correctly.	Inspect lamp installation.
	Lamp is defective.	Replace lamp.
	Lamp circuit breaker in building is tripped.	Reset the lamp circuit breaker.

PROBLEM	CAUSE	CORRECTIVE ACTION
Fume hood blower operates but lights will not operate	Lamp wiring is disconnected.	Inspect lamp wiring.
	Defective lamp ballasts.	Replace lamp ballasts.
Contaminants outside of fume hood	Improper user techniques for the fume hood.	See “Certifying the Hood” Chapter 3 and “Safety Precautions” Chapter 4 sections in the manual. (Ref. Appendix E)
	Restriction of the baffle air slots or – blockage of the exhaust outlet.	Remove baffles to ensure that all air slots, and the exhaust outlet are unobstructed.
	External factors are disrupting the fume hood airflow patterns or acting as a source of contamination.	See “Location Requirements” Chapter 2, “Certifying the Hood” Chapter 3, and “Safety Precautions” Chapter 4 sections of this manual. (Ref. Appendix E)
	Fume hood has improper face velocity.	Have fume hood re-certified and check remote blower exhaust system. Hood should have average face velocity of 60-100 fpm.
Vertical sash no longer operates smoothly	Cable is frayed or plastic protection is damaged.	Inspect cable and replace cable if worn or damaged immediately; otherwise injury could result.
	Pulley bearing is damaged.	Replace pulley, bearing or add grease.
	Cable or chain has slipped off the pulleys or sprockets.	Re-install, cable or chain must be replaced immediately if damaged.
	Weight has broken pulleys or sprockets.	Replace pulleys or sprockets.
Combination A-Style sash no longer operates smoothly	Horizontal glass panels have come off the tracks.	Re-install horizontal glass on tracks.

Chapter 8: Troubleshooting

PROBLEM	CAUSE	CORRECTIVE ACTION
	Vertical sash frame is distorted.	Place horizontal glass symmetrically and pull sash down to air foil. Straighten damaged frame.
	Cable is frayed or has slipped off the pulleys.	Re-install, cable must be replaced immediately if damaged.
Electrical duplex outlets no longer have power	Wires not connected or faulty duplex.	Check wire connection or replace duplex.
	Circuit breakers tripped in building electrical supply.	Reset circuit breakers.
Service valves no longer operate	Faulty building supply.	Inspect building supply shut off valves and appropriate pressures below the recommended 40 PSI.
	Valve no longer operates.	Replace valve and check for leaks.
	Supply line or outlet line has leaks.	Inspect line for leaks and fix any leaking plumbing connections.

APPENDIX A

PROTECTOR XL COMPONENTS

Illustration A-1 indicate the location of the following service parts:

Protector XL Replacement Parts

Item	Quantity	Part No.	Description
1A	1	9823700	Valve, Labconco (Water) 1/4" Compression Fitting
1B	1	9823701	Valve, Labconco (Water) 3/8" Compression Fitting
1C	1	9817000	Valve, Labconco 1/4" Compression Fitting (AIR, GAS, VAC, NIT, etc.)
1D	1	9817001	Valve, Labconco 3/8" Compression Fitting (AIR, GAS, VAC, NIT, etc.)
1E	1	9823702	Valve, Labconco Deionized 1/4" Compression Fitting
1F	1	9823703	Valve, Labconco Deionized 3/8" Compression Fitting
1G	1	9818000	Nut, Valve Mounting (Labconco)
2A	1	9826800	WaterSaver Valve/Gooseneck – GRN
2B	1	9826801	WaterSaver Valve/Connector (VAC) – YEL
2C	1	9826802	WaterSaver Valve/Connector (AIR) – ORG
2D	1	9826803	WaterSaver Valve/Connector (GAS) – BLU
2E	1	9826805	WaterSaver Valve/Connector (HOT WATER) – RED
2F	1	9826806	WaterSaver Valve/Connector (CW) – GRN
2G	1	9826807	WaterSaver Valve/Connector (STEAM) – BLK
2H	1	9826808	WaterSaver Valve/Connector (NITROGEN) – BRN
2I	1	9826809	WaterSaver Valve/Connector (OXYGEN) – LIGHT GREEN
2J	1	9826810	Swivel Gooseneck only – GRN
2K	1	9826812	Swivel Gooseneck only – WHITE
3	1	9818700 thru 08	Knobs (GRAY, GRN, BLU, ORG, YEL, RED, WHT, BLK, BRN)
4A	1	9818800	Hose Barb, GRAY (NEUTRAL OR ARGON) – NOT SHOWN
4B	1	9818801	Hose Barb, GREEN (COLD WATER) – NOT SHOWN
4C	1	9818802	Hose Barb, BLUE (GAS) – NOT SHOWN
4D	1	9818803	Hose Barb, ORANGE (AIR) – NOT SHOWN
4E	1	9818804	Hose Barb, YELLOW (VACUUM) – NOT SHOWN
4F	1	9818805	Hose Barb, RED (HOT WATER) – NOT SHOWN
4G	1	9818806	Hose Barb, WHITE (DEIONIZED WATER) – NOT SHOWN
4H	1	9818807	Hose Barb, BLACK (NEUTRAL OR STEAM) – NOT SHOWN
4I	1	9818808	Hose Barb, BROWN (NITROGEN) – NOT SHOWN
4J	1	9819000	Nut, Hose Barb – NOT SHOWN
5	1	9825500	Label, Knob (contains all the labels)
6	1	9818900	Lens, Knob
7A	1	9947100, 01, 02	115V Duplex Receptacle (GRAY) Right, Left 4' - 6', Left 8' w/ wires
7B	1	9818200	Cover Plate 115V Duplex
7C	1	9947103, 04, 05	115V GFCI Duplex Receptacle (GRAY) Right 4' - 6', Left 8' w/ wires
7D	1	9818100	Cover Plate, 115V GFCI
7E	1	9818300	Cover Plate, Blank
7F	1	9851203	115V Duplex Receptacle, Left 10' - 12' (GRAY)
7G	1	9851303	115V Duplex Receptacle, Left 16' (GRAY)
7H	1	9851603	115V GFCI Duplex Receptacle, Left 10' - 12' (GRAY)
7I	1	9851703	115V GFCI Duplex Receptacle, Left 16' (GRAY)

Appendix A: Protector XL Components

Item	Quantity	Part No.	Description
8A	1	1297503	LED, Direct Wired T8 x 3' – newest models use on 4', 8' & 16' Hoods – 12 watts each
8B	1	1297504	LED, Direct Wired T8 x 4' – newest models use on 5', 6', 10' & 12' Hoods – 22 watts each
8C	1	9945300	LED Light Fixture, 36" – use on 4' Hoods
8D	1	9945301	LED Light Fixture, 48" – use on 5', 6' and 7' Hoods
8E	1	9945302	LED Light Fixture, 36" Combo – use on 8' Hoods
8F	1	9945100, 01, 02	LED Light Fixture, Combo, use on 10', 12' and 16' Hoods
10A	1	1302301	Switch, Rocker
10B	1	1327500	Switch, Plug (Fills cutout when switch is not used)
11	1	9818400	Access Cover
12A	1	9409801	Side Panel, 30" internal deep hoods, upper
12B	1	9409802	Side Panel, 36" internal deep hoods, upper
12C	1	9409803	Side Panel, 48" internal deep hoods, upper
12D	4	1885308	Screw, Machine #10-24 x .50 Phillips
13A	1	9477400	4' Vertical Sash, Lower
13B	1	9441301	4' Vertical Sash, Upper
13C	1	9477401	5' Vertical Sash, Lower
13D	1	9441302	5' Vertical Sash, Upper
13E	1	9477402	6' Vertical Sash, Lower
13F	1	9441303	6' Vertical Sash, Upper
13G	1	9477403	8' Vertical Sash, Lower
13H	1	9441305	8' Vertical Sash, Upper
13I	4	9716604	8' Horizontal Sash
13J	4	9716605	10' Horizontal Sash
13K	4	9716606	12' Horizontal Sash
13L	4	9716607	16' Horizontal Sash
14	2	1861400 or 1850000	Pulley, Front, 1-3/16 Dia. or 2" Dia. Nylon
15A	2	4949902	Cable, Sash 130" – Upper Sash – NOT SHOWN
15B	2	4949904	Cable, Sash 190" – Lower Sash
15C	2	9414013	Cable Replacement Kit, 30" Interior Walk-In
15D	2	9414014	Cable Replacement Kit, 36" Interior Walk-In
15E	2	9414015	Cable Replacement Kit, 48" Interior Walk-In
15F	1	9545800	Weight Support Bracket Kit
16	2	9709300	Plastic Pulley, (Rear 2")
17	2	9713300	Bumper, Rubber (lower sash bumper)
18	4	1934100	Bronze Bearing, Flanged Rear
19	4	1920100	Clamp, Cable Replacement – NOT SHOWN
20	2	9935800	Threaded Connector (to attach weight to cable) – NOT SHOWN
21	1	9946300 or 01	Wiring Harness, Main
22A	1	9409600, 01, 02, 03, 04, 05	Front Panel, 3', 4', 5', 6', 7', 8'
22B	1	9499500, 01, 02, 03, 04, 05	Front Panel, 10' L, 10' R, 12' L, 12' R, 16' L, and 16' R
23A	1	9543900	Side Panel, 30" internal deep, lower
23B	1	9543902	Side Panel, 36" internal deep, lower
23C	1	9543904	Side Panel, 48" internal deep, lower

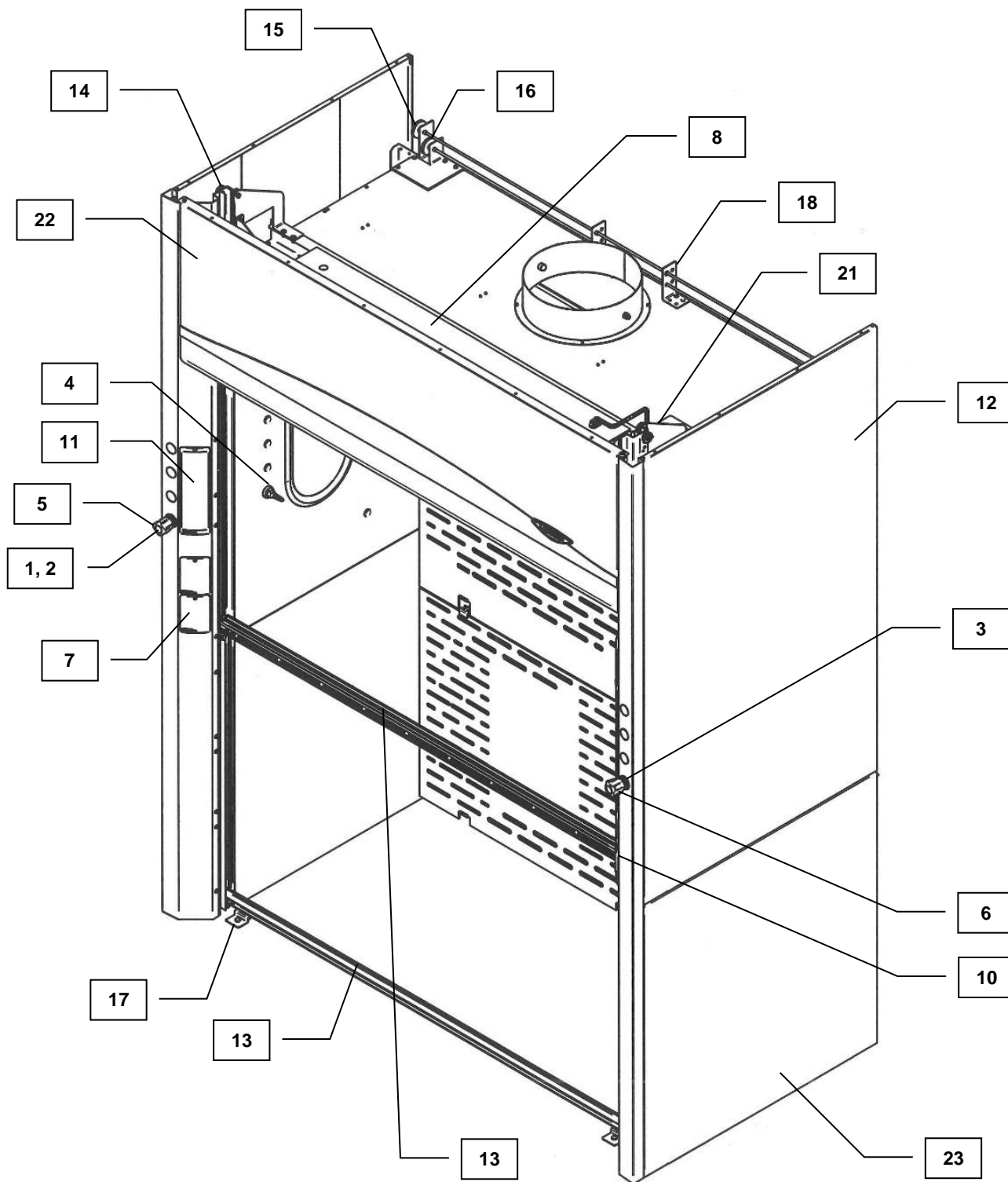
Figure A-1
Vertical Replacement Parts

Appendix A: Protector XL Components

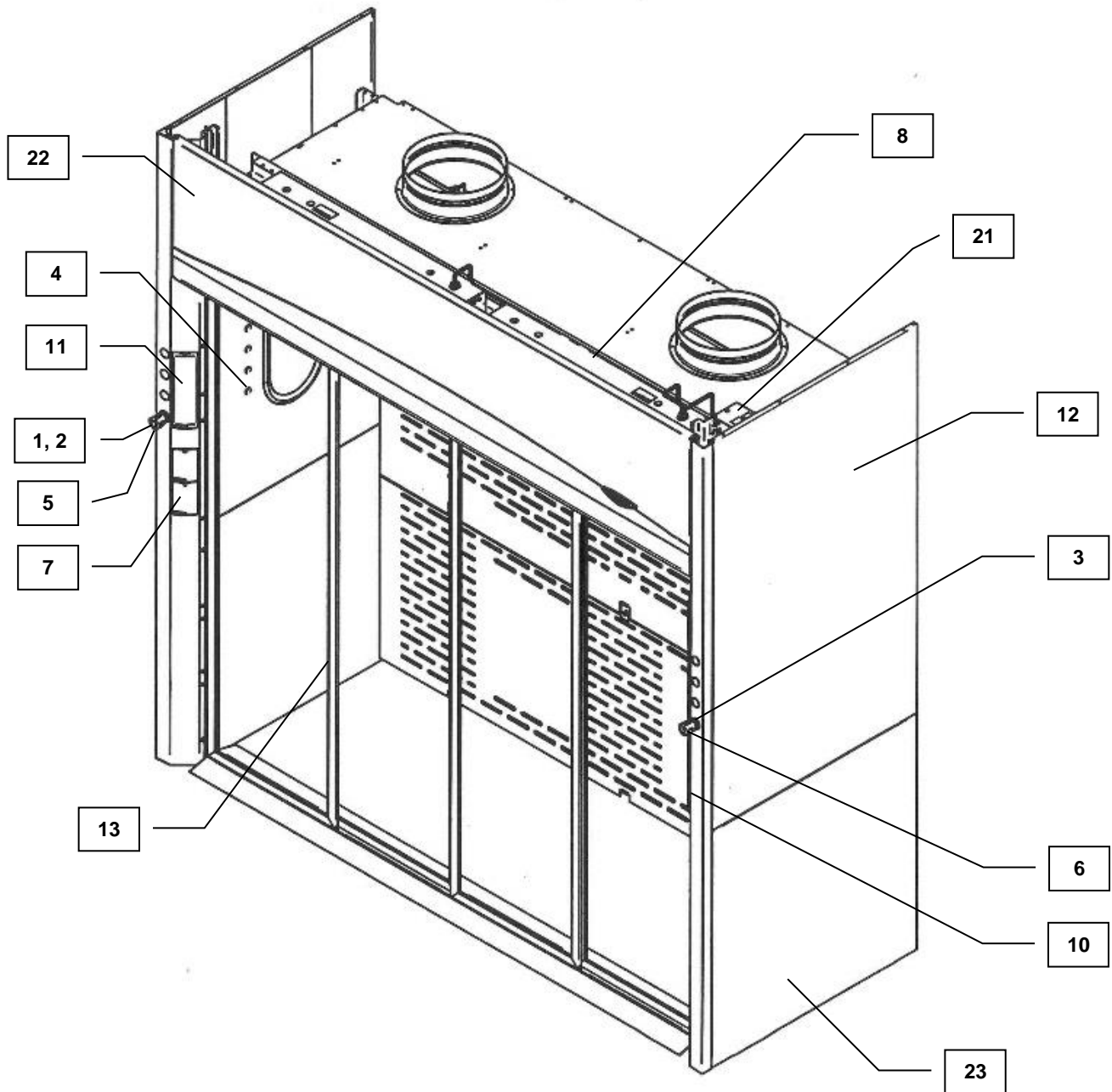
Figure A-2
Horizontal Replacement Parts

APPENDIX B

PROTECTOR XL DIMENSIONS

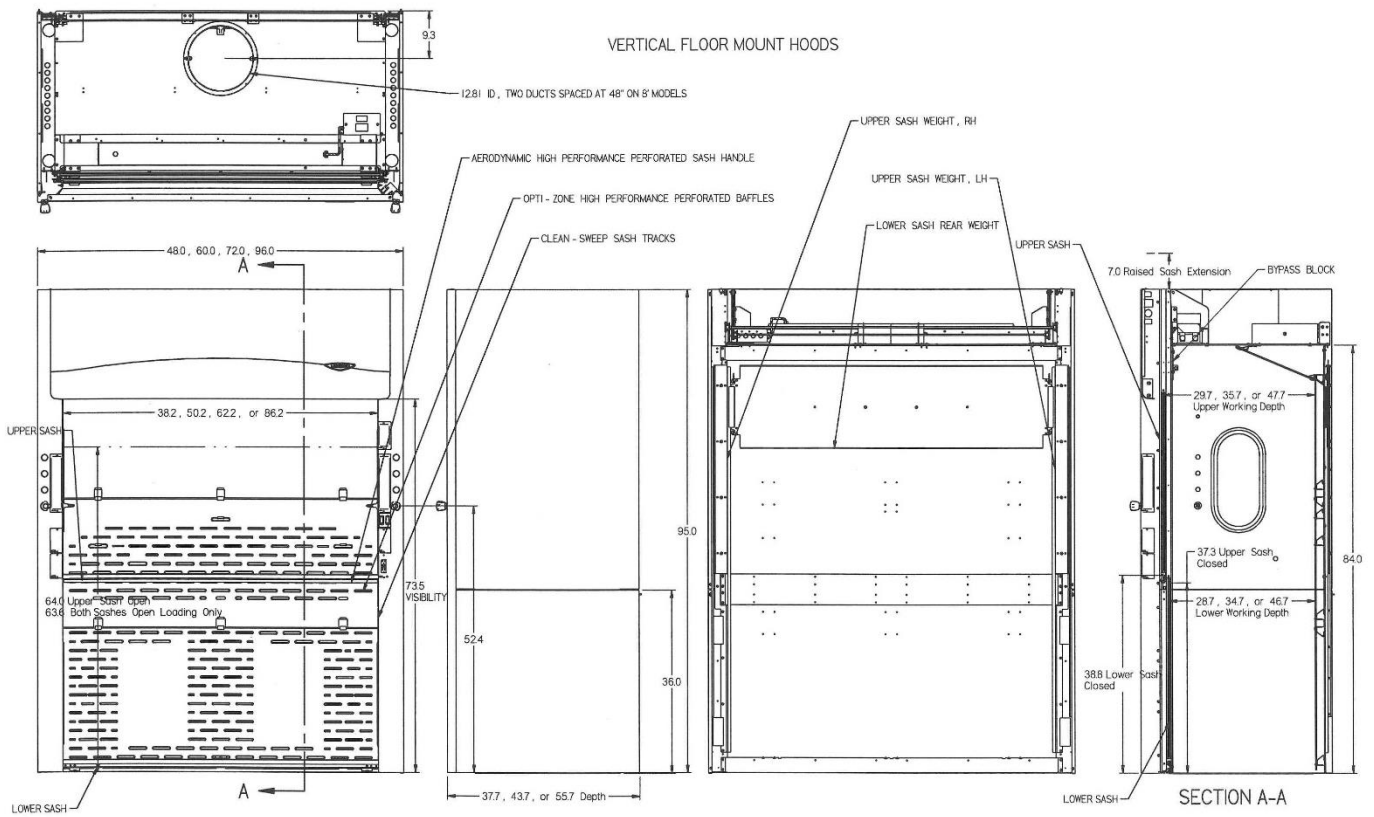
Figure B-1
Vertical Floor-Mount Dimensions

Appendix B: Protector XL Dimensions

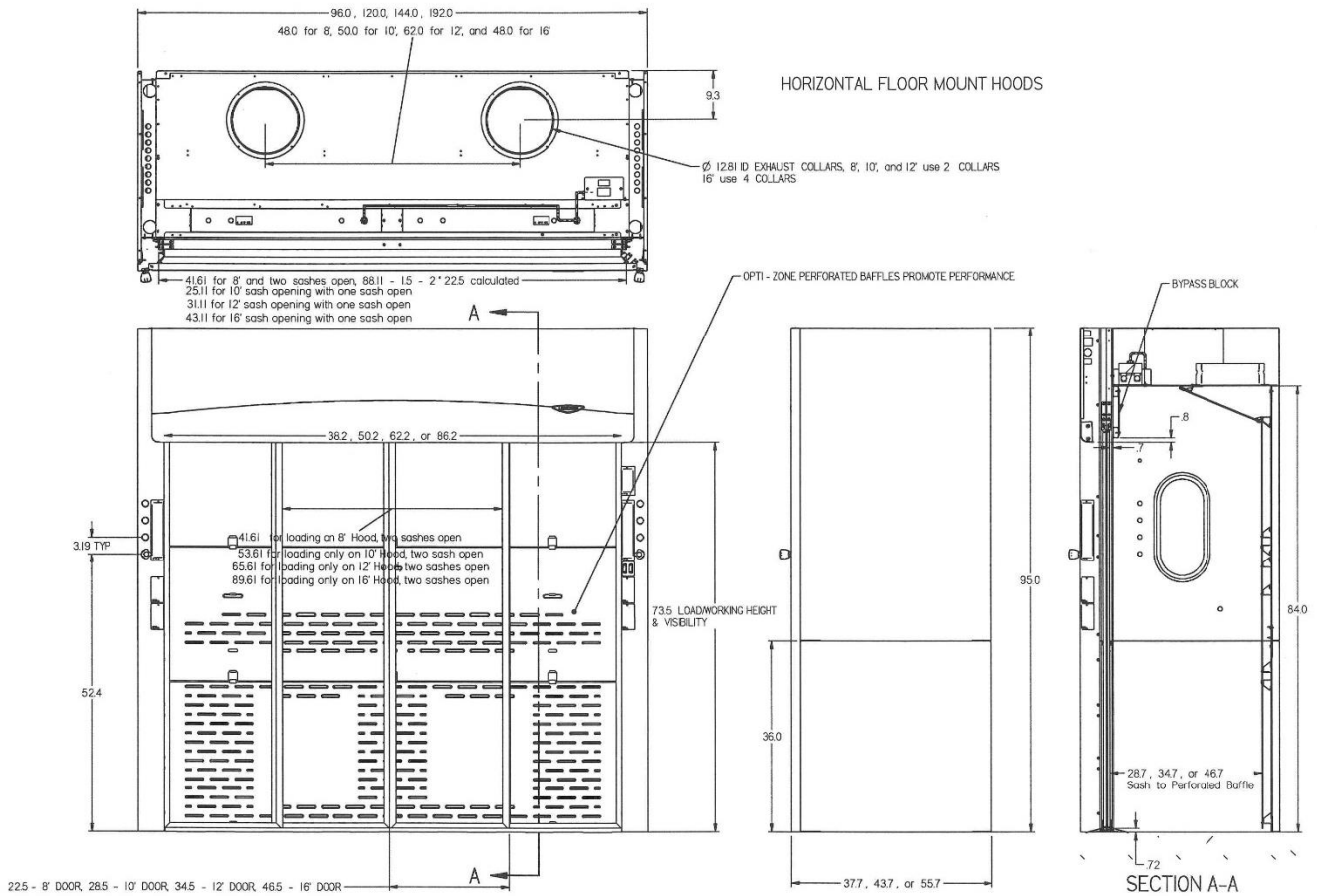
Figure B-2
Horizontal Floor-Mount Dimensions

APPENDIX C

PROTECTOR XL

SPECIFICATIONS


Environmental Conditions

- Indoor use only.
- Maximum altitude: 10,000 feet (3,048 meters).
- Ambient temperature range: 41° to 104°F (5° to 40°C).
- Maximum relative humidity: 80% for temperatures up to 88°F (31°C), decreasing linearly to 50% relative humidity at 104°F (40°C).
- Main supply voltage fluctuations not to exceed $\pm 10\%$ of the nominal voltage.
- Transient over-voltages according to Installation Categories II (Over-voltage Categories per IEC 1010). Temporary voltage spikes on the AC input line that may be as high as 1500V for 115V models and 2500V for 230V models are allowed.
- Used in an environment of Pollution degrees 2 (i.e., where normally only non-conductive atmospheres are present). Occasionally, however, a temporary conductivity caused by condensation must be expected, in accordance with IEC 664.
- Electrical Ratings


<u>Volts AC</u>	<u>Phase</u>	<u>Cycle</u>	<u>AMP*</u>
115	1	50/60	10
115	1	50	10
115	1	60	10
230	1	50/60	5
230	1	50	5
230	1	60	5

*Does not include current rating of receptacles.

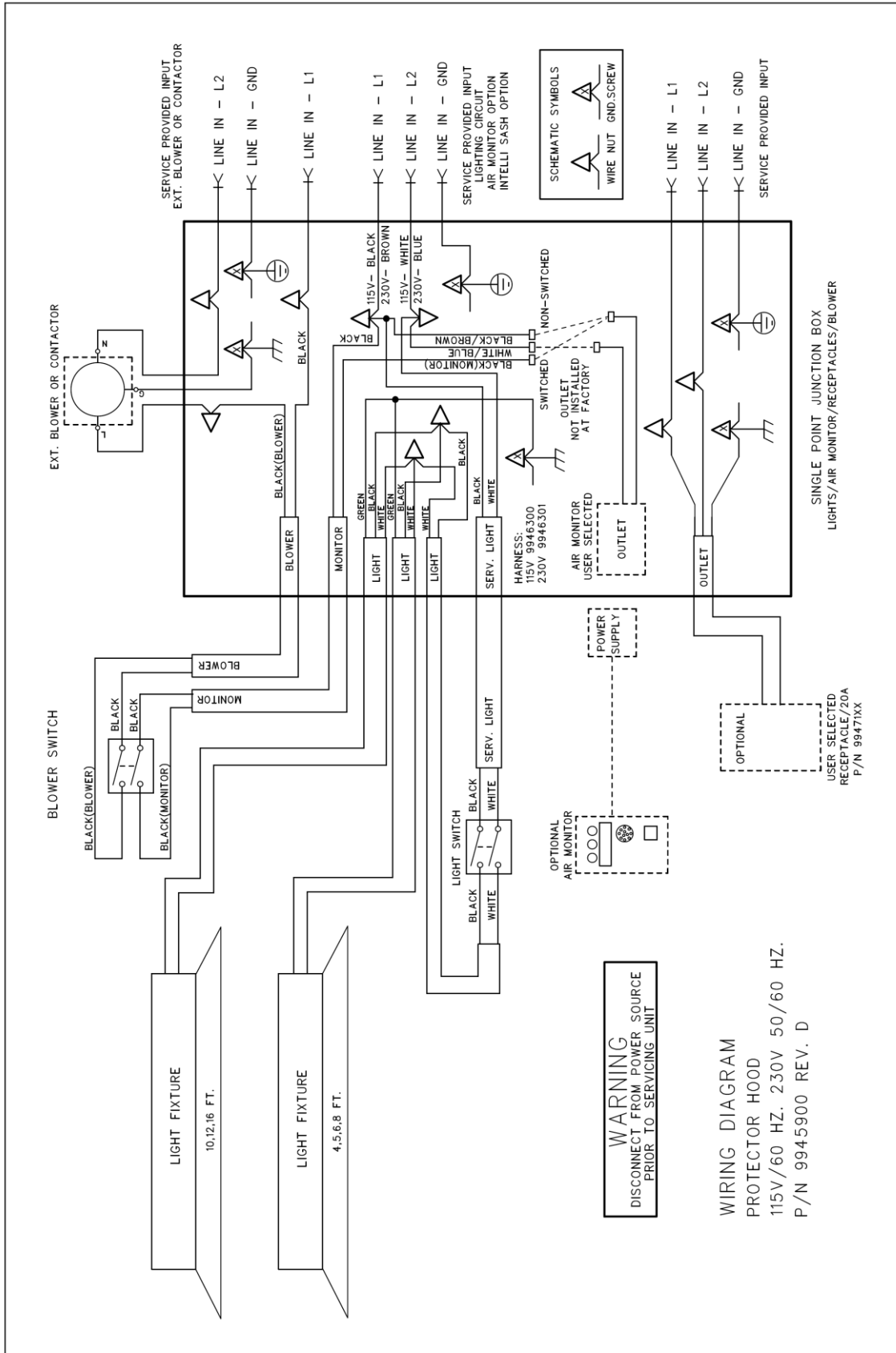Appendix C: Protector XL Specifications

Appendix C: Protector XL Specifications

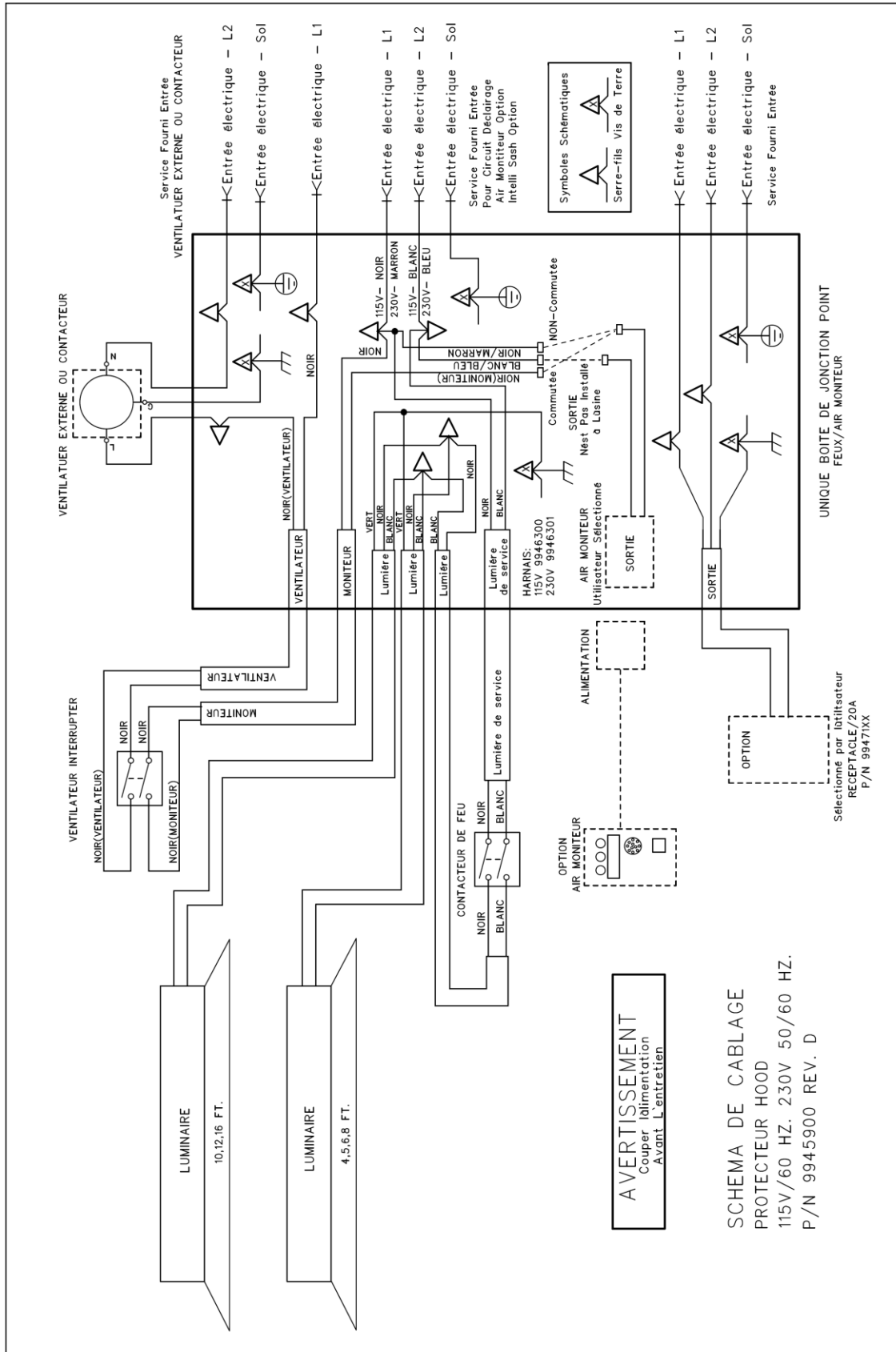APPENDIX D

SERIAL NUMBER TAG

DESCRIPTION

Serial tag includes standard information with the following changes:

The portion of the tag for electrical information is labeled with the following units: Volts AC, Phase, Cycle, AMP.

Each hood will include one of the following:

<u>Volts AC</u>	<u>Phase</u>	<u>Cycle</u>	<u>AM</u>  <u>X</u>
115	1	50/60	10X
115	1	50	10X
115	1	60	10X
230	1	50/60	5X
230	1	50	5X
230	1	60	5X

Where  may be a separate label indicating: Caution, See Manual.

The X above will be one of several possible alpha characters as defined in the manual as follows:

- X: No additional duplexes on hood, noted amperage applies.
- B: This fume hood includes **one** additional electrical receptacle, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.
- C: This fume hood includes **two** additional electrical receptacles, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.

- D: This fume hood includes **three** additional electrical receptacles, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.
- E: This fume hood includes **four** additional electrical receptacles, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.
- F: This fume hood includes **one** additional electrical receptacle, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- G: This fume hood includes **two** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- H: This fume hood includes **three** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- J: This fume hood includes **four** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- K: This fume hood includes **one** additional electrical receptacle, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- L: This fume hood includes **two** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.

Appendix D: Serial Number Tag Description

- M: This fume hood includes **three** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- N: This fume hood includes **four** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.

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APPENDIX E

REFERENCES

Many excellent reference texts and booklets are currently available. The following is a brief listing:

Laboratory Ventilation Standards

Federal Register 29 CFR Part 1910

Non-mandatory recommendations from "Prudent Practices".

- Fume hoods should have a continuous monitoring device
- Face velocities should be between 60-100 linear feet per minute (lfpm)
- Average 2.5 linear feet of hood space per person

Occupational Health and Safety

U.S. Department of Labor

200 Constitution Avenue N.W.

Washington, DC 20210

(202) 523-1452

Industrial Ventilation-ACGIH

- Fume hood face velocities between 60-100 lfpm
- Maximum of 125 lfpm for radioisotope hoods
- Duct velocities of 1000-2000 fpm for vapors, gasses and smoke
- Stack discharge height 1.3-2.0 x building height
- Well designed fume hood containment loss, <0.10 ppm

Industrial Ventilation, A Manual of Recommended Practice.

24th Edition, 2001

American Conference of Governmental Industrial Hygienists

1330 Kemper Meadow drive

Cincinnati, OH 45240-1634

(513) 742-2020

ASHRAE 110-1995 Method of Testing Performance of Fume Hoods

Evaluates fume hood's containment characteristics

- Three part test: Smoke generation, Face velocity profile, Tracer gas release @ 4 liters per minute
- Rated As Manufactured (AM), As Installed (AI) and As Used (AU)

American Society of Heating, Refrigerating, and Air Conditioning Engineers

1791 Tullie Circle N.E.

Atlanta, GA 30329

(404) 636-8400

ANSI Z9.5-2011 Laboratory Standard

Covers entire laboratory ventilation system.

- New and remodeled hoods shall have a monitoring device
- Ductless hoods should only be used with non-hazardous materials

American Industrial Hygiene Association

2700 Prosperity Avenue, Suite 250

Fairfax, VA 22031

(703) 849-8888

SEFA 1-2002

- Fume hood face velocities based on toxicity levels of chemicals
 - Class A – 125 to 150 fpm
 - Class B – 80 to 100 fpm
 - Class C – 75 to 80 fpm
- Test method – face velocity profile and smoke generation

Scientific Equipment & Furniture Association

1028 Duchess Drive

McLean, VA 22102

(703) 538-6007

NFPA 45 – 2002 Fire Protection for Laboratories Using Chemicals

- Laboratory hoods should not be relied on for explosion protection
- Exhaust air from fume hoods should not be recirculated
- Services should be external to the hood
- Canopy hoods only for non-hazardous applications
- Materials of construction should have flame spread of 25 or less
- 80 to 120 fpm to prevent escape

NFPA 30 – 2000 Flammable and Combustible Liquids Code

- Approved cabinets may be metal or wood
- Vent location on cabinets are required
- Venting of cabinets not a requirement

National Fire Protection Association

1 Batterymarch Park

P.O. Box 9101

Quincy, MA 02269-9101

(800) 344-3555

General References

American Conference of Governmental Industrial Hygienists. *Industrial Ventilation, A Manual of Recommended Practice*, Cincinnati, OH

ASHRAE Standard Committee. *ASHRAE Standard* Atlanta: ASHRAE Publications Sales Department, 1995

British Standards Institution, *Laboratory Fume Cupboards*. Parts 1, 2 and 3, London: 1990

Department of Labor, Occupational Safety and Health Administration, *29 CFR Part 1910, Occupational Exposures to Hazardous Chemicals in Laboratories, Final Rule*. Vol. 55, No. 21. Washington D.C.:1990

Appendix E: References

DiBerardinis, L. et al. *Guides for Laboratory Design, Health and Safety Considerations*. Wiley & Sons, 1987

McDermott, Henry, *Handbook of Ventilation for Contaminant Control, 2nd Edition*. Butterworth Publishers, 1985.

Miller, Brinton M. et al. *Laboratory Safety: Principles and Practices*. American Society for Microbiology, Washington, D.C.: 1986

NIH Guidelines for the Laboratory Use of Chemical Carcinogens. NIH Publication No. 81-2385.

Rayburn, Stephen R. *The Foundation of Laboratory Safety, A Guide for the Biomedical Laboratory*. Springer-Verlag, New York: 1990

Sax, N. Irving and Lewis, JR., Richard J. *Rapid Guide to Hazardous Chemicals in the Workplace*. Van Nostrand Reinhold, 1987.

Schilt, Alfred A. *Perchloric Acid and Perchlorates*. The G. Frederick Smith Chemical Company, Columbus, OH: 1979.

Steere, Norman. *CRC Handbook of Laboratory Safety, 2nd Edition*. CRC Press, 1971.

APPENDIX F

DISASSEMBLY AND REASSEMBLY INSTRUCTIONS

Required Tools

The following tools will be required for disassembly and reassembly:

Cordless Drills and Phillips bits.

1/4", 5/16", 3/8", 7/16" Sockets	Small Screwdriver
Phillips Screwdriver	Right Angle Screwdriver
1/4", 5/16" Box Ratchet	1/4" Extension
5/16", 3/8", 7/16" Combination Wrench	1/4" Ratchet
Small Right Angle Needle Nose Pliers	Nut Driver Set

Decorative Panel Removal

1. Remove side panels by removing #10-24 screws and lifting away.
2. Remove front panels by lifting up and away.

Light Removal and Blower Wiring

1. On the top of the hood, disconnect the Molex connector and remove the green ground wire. All the removable wires are part of the wiring harness to the corner post switches. Be sure to label all wires and refer to the wiring diagrams for proper reassembly.

Header Removal

1. The header is fastened to the corner posts by four #12 screws. Simply reach behind the corner posts from the side and remove the screws while another person supports the header.

Appendix F: Disassembly & Reassembly Instructions

Sash Removal

There is one sash to be removed on the 4', 5', and 6' hoods. Please note that there are two sashes that need to be removed for the 8', 10', and 12' hoods and four sashes on the 16' hood. Please repeat the procedure for each sash.

1. For 8' through 16', remove the bumper from the middle sash slide by removing the screw. Be careful not to misplace the loose parts. This part is a crucial component to the operation of the sash.
2. For 8' through 16', remove the middle sash slide assemblies by sliding them up and out.
3. Remove the sash end stops by removing the hardware.
4. Remove end cable and pulleys from sides.
5. Now the glass sash assembly can be removed by sliding it up and out.
6. Remove sash weights. The weights were shipped attached to the back.

Top Assembly Removal

Please note that the 4' through 8' hoods are designed with one continuous top liner. Therefore, the top support is not required for the 4' through 8' hoods, but is used on the 10', 12', and 16' hoods.

1. Remove the top support by removing the 1/4 –20 hardware and screws.
2. Remove the anti-rack shaft assemblies by removing the screws.
3. Then remove the top liner assemblies by removing the screws.

Front Truss Removal (if applicable on larger hoods 8' and up)

1. Remove the front truss support from the side frames by removing the 1/4 –20 x 1.5" hex head screws, washers, lockwashers, and nuts. Simply reach behind the side frames while somebody supports the front truss and remove it.

Airfoil Removal

1. Remove lower corner post screw located under the airfoil.
2. Remove the airfoil by removing the rear screws located on top of the airfoil on each side.

Rear Frame Removal

1. Remove the rear frames from the side frames by removing screws from each side frame.
2. If desired, you can remove the right rear frame from the left rear frame by removing the 1/4-20 hardware, angle support, and screws. This step is not always necessary and depends on the installation and size of the hoods.

Transport and Reassembly

The various subassemblies are now ready to be transported and reassembled in the opposite way the hood was disassembled. All hoods will require a minimum of five people to lift the hood, two people on each side and one in front (for balance only). A hydraulic lift table should be used as an aid in lifting. The 16' hood should be assembled where placed. If in doubt about lifting the hood, always reassemble the hood where it will be placed.

Reinstallation

Make sure the anti-rack shaft does not slide from left to right as this will be noisy and not work effectively.

Sash Weight Reassembly

After the hood is assembled, the sash weights are easily installed before the hood is placed against a rear wall. **Please keep this in mind and install the sash weights now.**

Baffles Installation

Install the upper baffle, middle baffle, and lower baffle. Be sure the baffles are resting in the proper baffle mount supports.

Hood Replacement and Service Installation

The hood is now ready for placement and the HVAC, plumbing and electrical service can be installed.

APPENDIX G

VAV FLOOR-MOUNTED

OPEN AREA GUIDELINES

When integrating a Labconco floor-mounted fume hood into a variable air volume (VAV) mechanical system, properly assessing the area opening of the fume hood is critical to ensure correct controller calibration for maintaining face velocity across all sash positions. Depending on your brand of VAV controller, you may be prompted to enter different information to arrive at this correct area. To simplify this requirement, Labconco offers the following step-by-step procedure to ensure you have the correct information.

1. Locate the serial tag for your floor-mounted fume hood; usually on the lower right of the sash glass.
2. Move to the section of the document below specific to your sash type and model number.

Model numbers beginning 112

Move to the section below specific to your sash type

Vertical rising sashes (two sashes that move up and down)

1. Determine the free opening area
 - a. Width
 - i. Measure the interior width of the fume hood just behind the plane of the sash glass and enter that here:

Width Opening _____ inches = A

- b. Height
 - i. Measure the vertical distance between the top of the closed lower sash and the bottom of the sash handle on the upper sash, and enter that here:

Vertical Opening _____ inches = B

Appendix G: VAV Floor-Mounted Open Area Guidelines

- c. Free opening area
- i. Multiply the width opening times the vertical opening (A x B), and enter that here:

Free opening area = _____ square inches = C (sq in)

- ii. Divide this by 144 to arrive at square feet, and enter that here:

Free opening area = _____ square feet = C (sq ft)

2. Determining bypass or “leakage” area

- a. This is the opening width (A) times 2.2 inches (A x 2.2 = bypass area).

Bypass area = _____ square inches = D (sq in)

- b. Divide this by 144 to arrive at square feet

Bypass area = _____ square feet = D (sq ft)

3. Determining Total Effective Opening Area

- a. This is the Free Opening Area plus the Bypass Opening Area (C + D = Total Effective Opening Area)

Total Effective Area = _____ square inches

Total Effective Area = _____ square feet

Model numbers beginning 112 (continued)

Horizontal sashes (Sashes that move left and right)

1. Determine the free opening area

- a. Width

- i. Measure the opening between sashes and enter that here:

Width Opening _____ inches = A

- b. Height

- i. Measure the vertical distance between the top of the sash track and bottom of the header panel, and enter that here:

Vertical Opening _____ inches = B

Appendix G: VAV Floor-Mounted Open Area Guidelines

- c. Free opening area
- i. Multiply the width opening times the vertical opening (A x B), and enter that here:
Free opening area = _____ square inches = C (sq in)
 - ii. Divide this by 144 to arrive at square feet, and enter that here:
Free opening area = _____ square feet = C (sq ft)

2. Determine bypass or “leakage” area (D) from the table below

Nominal exterior width (ft)	Bypass or "leakage" area (sq ft)	Bypass or "leakage" area (sq in)
8	0.47	68
10	0.96	138
12	1.07	154
16	1.35	194

3. Determining Total Effective Opening Area
- a. This is the Free Opening Area plus the Bypass Opening Area (C + D = Total Effective Opening Area)
Total Effective Area = _____ square inches
Total Effective Area = _____ square feet